




Product Information

SIGMA QUALITY CONTROL TEST PROCEDURE

Enzymatic Assay of CHYMOTRYPSIN¹ (EC 3.4.21.1)

PRINCIPLE:

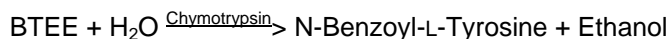

Abbreviation used:

BTEE = N-Benzoyl-L-Tyrosine Ethyl Ester

CONDITIONS: T = 25°C, pH = 7.8, $A_{256\text{nm}}$, Light path = 1 cm

METHOD: Continuous Spectrophotometric Rate Determination

REAGENTS:

- A. 80 mM Tris HCl Buffer, pH 7.8 at 25°C
(Prepare 100 ml in deionized water using Trizma Base, Sigma Prod. No. T-1503. Adjust to pH 7.8 at 25°C with 1 M HCl.)
- B. 1.18 mM N-Benzoyl-L-Tyrosine Ethyl Ester Solution (BTEE)
(Prepare 50 ml using N-Benzoyl-L-Tyrosine Ethyl Ester, Sigma Prod. No. B-6125, by initially dissolving in 31.7 ml of Methanol, Absolute, Sigma Stock No. 17-5. Add enough deionized water to make the final volume 50 ml.)
- C. 2 M Calcium Chloride Solution (CaCl_2)
(Prepare 5 ml in deionized water using Calcium Chloride, Dihydrate, Sigma Prod. No. C-3881.)
- D. 1 mM Hydrochloric Acid Solution (HCl)
(Prepare 50 ml in deionized water using Hydrochloric Acid, Sigma Prod. No. H-7020.)
- E. Chymotrypsin Enzyme Solution
(Immediately before use, prepare a solution containing 2 - 5 units/ml of Chymotrypsin in cold Reagent D.)

Enzymatic Assay of CHYMOTRYPSIN¹
(EC 3.4.21.1)

PROCEDURE:

Pipette (in milliliters) the following reagents into suitable quartz cuvettes:

	<u>Test</u>	<u>Blank</u>
Reagent A (Buffer)	1.42	1.42
Reagent B (BTEE)	1.40	1.40
Reagent C (CaCl ₂)	0.08	0.08

Mix by inversion and equilibrate to 25°C. Monitor the A_{256nm} until constant, using a suitably thermostatted spectrophotometer. Then add:

Reagent D (HCl)	-----	0.10
Reagent E (Enzyme Solution)	0.10	-----

Immediately mix by inversion and record the increase in A_{256nm} for approximately 5 minutes. Obtain the $\Delta A_{256nm}/\text{minute}$ using the maximum linear rate for both the Test and Blank.

CALCULATIONS:

$$\text{Units/ml enzyme} = \frac{(\Delta A_{256nm}/\text{min Test} - \Delta A_{256nm}/\text{min Blank})(3)(df)}{(0.964) (0.1)}$$

3 = Volume (in milliliters) of assay

df = Dilution factor

0.964 = Millimolar extinction coefficient of N-Benzoyl-L-Tyrosine at 256 nm.

0.1 = Volume (in milliliters) of enzyme used

$$\text{Units/mg solid} = \frac{\text{units/ml enzyme}}{\text{mg solid/ml enzyme}}$$

$$\text{Units/mg protein} = \frac{\text{units/ml enzyme}}{\text{mg protein/ml enzyme}}$$

UNIT DEFINITION:

One unit will hydrolyze 1.0 μmole of BTEE per minute at pH 7.8 at 25°C.

Enzymatic Assay of CHYMOTRYPSIN¹
(EC 3.4.21.1)

FINAL ASSAY CONCENTRATION:

In a 3.00 ml reaction mix, the final concentrations are 38 mM Tris, 0.55 mM N-benzoyl-L-tyrosine ethyl ester, 30% (v/v) methanol, 53 mM calcium chloride, 0.03 mM hydrochloric acid, and 0.2 - 0.5 unit chymotrypsin.

REFERENCES:

Wirnt, R. (1974) in *Methods of Enzymatic Analysis*, 2nd ed., Volume II, pp 1009-1012, Academic Press Inc., New York, NY

NOTES:

1. This assay procedure is not to be used to assay Chymotrypsin, Insoluble, Sigma Prod. Nos. C-9134 and C-7260, and Chymotrypsin, Acrylic Beads, Sigma Prod. No. C-5407.
2. This assay is based on the cited reference.
3. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that Sigma products conform to the information in Sigma publications. Purchaser must determine the suitability of the information and products for its particular use. Upon purchase of Sigma products, see reverse side of invoice or packing slip for additional terms and conditions of sale.