

Product Information

PERIODIC ACID-SCHIFF'S REAGENT

LEICA AutoStainer XL Stainer

(PAS-Light Green Template)
 Procedure 395

Station 1	Station 2	Station 3	Station 4	Station 5	Station 6	Station 7	W1	W2	W3	W4	W5	Oven
Xylene 1 min, 30 sec	Xylene 1 min, 30 sec	Xylene 1 min, 30 sec	100% Alc 1 min, 30 sec	100% Alc 1 min, 30 sec	95% Alc 30 sec	80% Alc 30 sec	Wash 5 sec	Wash 5 min	Wash 5 min	Wash 2 min	Wash 3 min, 30 sec	
End	Station 18 Xylene 2 min	Station 17 Xylene 2 min	Station 16 100% Alc 1 min	Station 15 100% Alc 20 sec	Station 14 100% Alc 10 sec	Station 13 95% Alc 5 sec	Station 12 Not Used	Station 11 Not Used	Station 10 Light Green 20 sec	Station 9 Schiff's Rgt 20 min	Station 8 Periodic Acid 5 min	Hold