



3050 Spruce Street
Saint Louis, Missouri 63103 USA
Telephone 800-325-5832 • (314) 771-5765
Fax (314) 286-7828
email: techserv@sia.com
sigma-aldrich.com

Product Information

Phosphate Buffered Saline, with BSA, pH 7.4

Product Number **P 3688**
Storage Temperature 2-8 °C

Product Description

This PBS-BSA formulation is recommended for use in immunohistology protocols.

Each pouch (3.5 in x 3.5 in) contains 8 g sodium chloride, 0.2 g potassium phosphate, monobasic, 1.5 g sodium phosphate, dibasic, 0.2 g potassium chloride, and 10.0 g BSA. There is 19.55 g/pouch. The contents of one pouch when dissolved in 1 liter of water, will yield 10 mM phosphate buffered saline (138 mM NaCl, 2.7 mM KCl), pH 7.4, at 25 °C with bovine serum albumin (1.0% w/v).

Precautions and Disclaimer

For Laboratory Use Only. Not for drug, household or other uses.

ALF/RXR 11/04

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.