


# OPTI-PRIME Priming Valve for Waters Pumps

## Product Specification

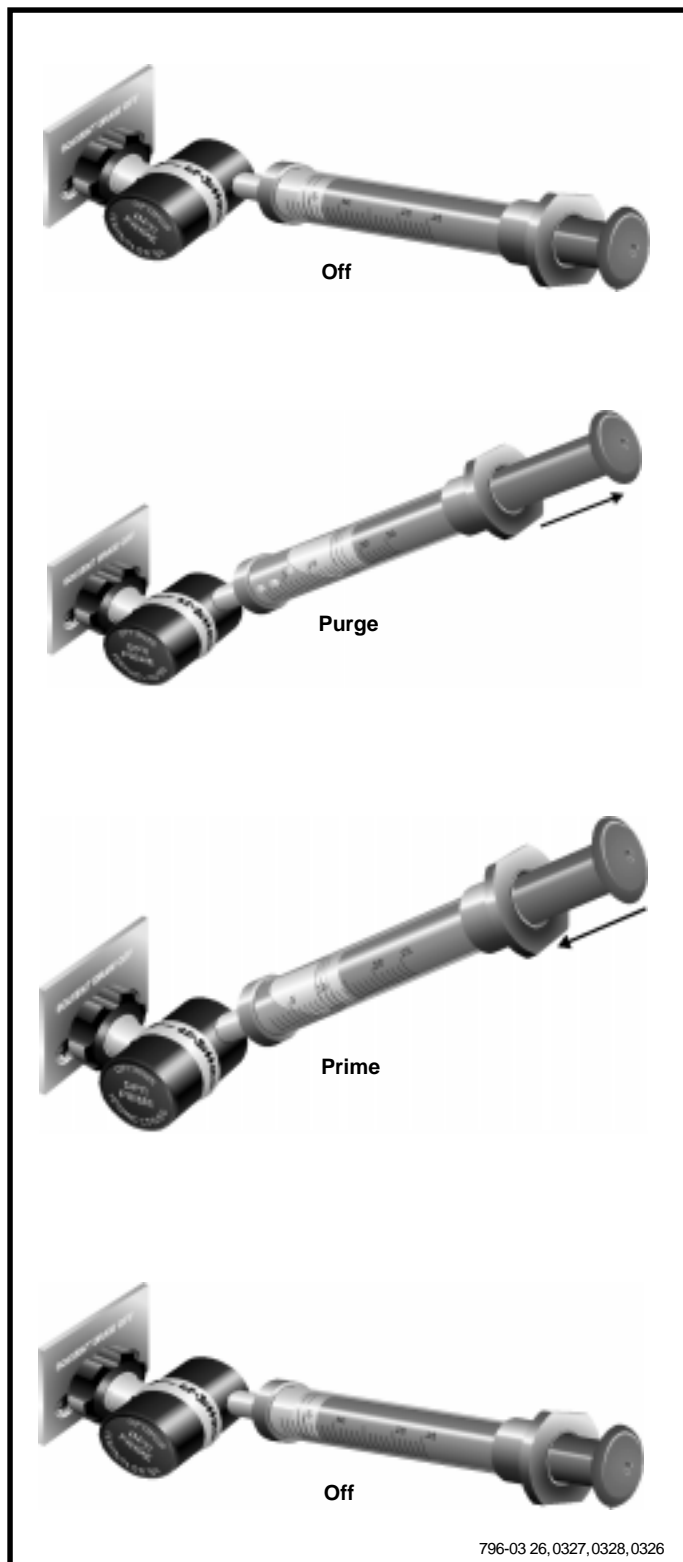

An OPTI-PRIME™ priming valve\* reduces the procedure for priming a Waters analytical pump to a simple task that takes only seconds. An easy, one-time installation eliminates wear and tear on the seal insert – there is no need to use the standard draw-off tube, nor to force a piece of Teflon® tubing over the stem.

### Benefits:

- *Single-handed operation*  
With the valve tightened in place and in the off (down) position, push the syringe into place on the valve. To open the valve, simply rotate the syringe upward.
- *No need to remove the syringe during priming*  
As you rotate the syringe upward to open the valve, air in the syringe will rise to the back of the barrel. This eliminates the need to remove the syringe to expel air.
- *No air is injected into the pump*  
Simply stop pushing on the syringe plunger before air reaches the valve. After priming the pump, rotate the syringe down to close the valve. If necessary, the syringe can be removed for refilling.

### Applications:

Use an OPTI-PRIME priming valve with Waters single solvent pumps M45, M45G, M501, M510, M590, M600E, M610, or with multiple solvent inlet pumps M6000 or M6000A.

### Construction:

Wetted surfaces are polyphenylene sulfide (PPS), polytetrafluoroethylene (PTFE), or 316 stainless steel, making the OPTI-PRIME valve inert to all commonly used HPLC solvents.

### Ordering Information:

Description	Cat. No.
OPTI-PRIME Priming Valve	59425-U
OPTI-PRIME Priming Valve with Optimize Safety Syringe	59426
Optimize Safety Syringe	59451

\*Patent pending.

### Trademarks

ChromFax — Supelco, Inc.

OPTI-PRIME — Optimize Technologies, Inc.

Teflon — E.I. du Pont de Nemours & Co., Inc.

## Free Technical Literature From Supelco

Supelco produces a wide variety of technical literature, all of which is free. You may request this literature several different ways:

- Use the business reply card located at the back of our catalog.
- Call our Ordering and Customer Service department (800-247-6628 or 814-359-3441).
- Use our ChromFax™ service.

### How to Use ChromFax

ChromFax is Supelco's automated digital data system, which allows you to instantly obtain technical literature any time of the day or night. Supelco technical bulletins, application notes, lab hints, and product specification sheets can be sent to your Fax machine.

1. To use ChromFax you must dial from a touch-tone phone. If you do not have a touch-tone phone, call our Ordering and Customer Service department and we will fax the publication to you.
2. Have the phone number of your Fax machine ready, and dial **800-652-8069 or 814-359-5748**. Remember you must have the *complete* Fax number (011 + country/city code + Fax number) when you are calling from outside the US or Canada, otherwise the document cannot be transmitted.

3. When ChromFax answers, follow the simple recorded instructions. Select the **ChromFax number** of the document you want. Be certain you order documents by their ChromFax numbers and NOT their literature numbers.
4. Enter your Fax number and hang up. The requested information will be sent to you within minutes. To report a problem with the ChromFax system contact our Technical Service department.

### Chemical Standards Data on ChromFax

ChromFax also allows you to receive analytical data for chemical standards. To receive this information, follow the ChromFax instructions outlined above. When ChromFax prompts you for the document number, enter the chemical standard's lot number, using digits only — do not enter alphabetical characters (e.g., lot number LA 1467: do not enter "LA").

### Technical Literature Available on ChromFax

The listing of documents available on ChromFax is updated every month. We encourage you to request a list of available ChromFax documents each time you use the ChromFax system.

---

For more information, or current prices, contact your nearest Supelco subsidiary listed below. To obtain further contact information, visit our website ([www.sigma-aldrich.com](http://www.sigma-aldrich.com)), see the Supelco catalog, or contact Supelco, Bellefonte, PA 16823-0048 USA.

**ARGENTINA** · Sigma-Aldrich de Argentina, S.A. · Buenos Aires 1119 **AUSTRALIA** · Sigma-Aldrich Pty. Ltd. · Castle Hill NSW 2154 **AUSTRIA** · Sigma-Aldrich Handels GmbH · A-1110 Wien  
**BELGIUM** · Sigma-Aldrich N.V./S.A. · B-2880 Bornem **BRAZIL** · Sigma-Aldrich Quimica Brasil Ltda. · 01239-010 São Paulo, SP **CANADA** · Sigma-Aldrich Canada, Ltd. · 2149 Winston Park Dr., Oakville, ON L6H 6J8  
**CZECH REPUBLIC** · Sigma-Aldrich s.r.o. · 186 00 Praha 8 **DENMARK** · Sigma-Aldrich Denmark A/S · DK-2665 Vallensbaek Strand **FINLAND** · Sigma-Aldrich Finland/YA-Kemia Oy · FIN-00700 Helsinki  
**FRANCE** · Sigma-Aldrich Chimie · 38297 Saint-Quentin-Fallavier Cedex **GERMANY** · Sigma-Aldrich Chemie GmbH · D-82041 Deisenhofen **GREECE** · Sigma-Aldrich (o.m.) Ltd. · Ilioupoli 16346, Athens  
**HUNGARY** · Sigma-Aldrich Kft. · H-1067 Budapest **INDIA** · Sigma-Aldrich Co. · Bangalore 560 048 **IRELAND** · Sigma-Aldrich Ireland Ltd. · Dublin 24 **ISRAEL** · Sigma Israel Chemicals Ltd. · Rehovot 76100  
**ITALY** · Sigma-Aldrich s.r.l. · 20151 Milano **JAPAN** · Sigma-Aldrich Japan K.K. · Chuo-ku, Tokyo 103 **KOREA** · Sigma-Aldrich Korea · Seoul **MALAYSIA** · Sigma-Aldrich (M) Sdn. Bhd. · 58200 Kuala Lumpur  
**MEXICO** · Sigma-Aldrich Quimica S.A. de C.V. · 50200 Toluca **NETHERLANDS** · Sigma-Aldrich Chemie BV · 3330 AA Zwijndrecht **NORWAY** · Sigma-Aldrich Norway · Torshov · N-0401 Oslo  
**POLAND** · Sigma-Aldrich Sp. z o.o. · 61-663 Poznań **PORTUGAL** · Sigma-Aldrich Quimica, S.A. · Sintra 2710 **RUSSIA** · Sigma-Aldrich Russia · Moscow 103062 **SINGAPORE** · Sigma-Aldrich Pte. Ltd.  
**SOUTH AFRICA** · Sigma-Aldrich (pty) Ltd. · Jet Park 1459 **SPAIN** · Sigma-Aldrich Quimica, S.A. · 28100 Alcobendas, Madrid **SWEDEN** · Sigma-Aldrich Sweden AB · 135 70 Stockholm  
**SWITZERLAND** · Supelco · CH-9471 Buchs **UNITED KINGDOM** · Sigma-Aldrich Company Ltd. · Poole, Dorset BH12 4QH  
**UNITED STATES** · Supelco · Supelco Park · Bellefonte, PA 16823-0048 · Phone 800-247-6628 or 814-359-3441 · Fax 800-447-3044 or 814-359-3044 · email: [supelco@sial.com](mailto:supelco@sial.com)

Supelco is a member of the Sigma-Aldrich family. Supelco products are sold through Sigma-Aldrich, Inc. Sigma-Aldrich warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product for a particular use. Additional terms and conditions may apply. Please see the reverse side of the invoice or packing slip.