

Product Information

Fetal Bovine Serum Sourced in USA heat inactivated

CATALOG NO. 12106

Description

Animal serum is commonly used to supplement basal media formulations for the optimal growth of many cell types *in vitro*.

Fetal Bovine Serum (FBS) is the most common serum used to supplement cell culture media due to its high nutritional content. Although it is relatively low in protein, FBS is effective in promoting and sustaining growth of vertebrate mammalian and insect cells.

Heat inactivation is commonly used to destroy complement. Heat inactivation is the method of choice of many investigators to insure that cells will not be lysed to antibody binding and complement activation. SAFC Biosciences' validated process consists of raising the temperature of the serum to 56 C for 30 minutes under controlled conditions.

Raw Serum Process

Animal blood is collected at United States Department of Agriculture (USDA) approved abattoirs located within the USA. Whole fetal blood is aseptically collected via cardiac puncture and allowed to clot under controlled conditions. After centrifugation, the serum is decanted from the clot. The raw serum is then pooled and immediately frozen. Raw serum is pre-tested by SAFC Biosciences according to established criteria, providing a complete awareness of all serum characteristics prior to filtration and further processing.

Filtration & Packaging

Frozen raw serum is thawed under controlled conditions and then processed through a series of membrane filters in descending pore size. Pooled FBS is filtered through three 0.1 µm filters. Integrity tests are conducted on the sterilizing filter pre- and post-filtration by bubble point and diffusive flow methods. The serum filtration process meets the sterility assurance level of 10⁻³ as verified by aseptic media fill validation.

Serum is dispensed under HEPA filtered, Class 100 conditions. Serum is packaged in sterilized, graduated plastic bottles and sealed with a tamper indicator. Bottles are identified with sequentially numbered labels and frozen at -10 to -40 C.

Traceability

The material used in this product is collected in the United States. Animals used for collection of serum are USDA inspected and acceptable for slaughter. A Certificate of Analysis indicating the country of origin is available for each lot of serum.

Precautions

This product is for further manufacturing use. THIS PRODUCT IS NOT INTENDED FOR HUMAN OR THERAPEUTIC USE. For stability and optimal performance, serum should be stored at -10 to -40 C and used prior to the labeled expiration date.

Use aseptic technique when handling serum. Refiltering sterile serum before or after being added to sterile medium is not recommended because the growth promoting capability may be reduced.

Serum may contain undetectable levels of adventitious agents. The risk of transmission of viruses can be minimized by the use of a validated virus inactivation process. Upon request, SAFC Biosciences will provide serum irradiated according to our validated SER-TAIN™ Gamma Irradiation process.

Storage

To effectively preserve the integrity of animal serum, it should be stored frozen and protected from light. For stability and optimal performance, serum should be stored at -10 to -40 C and used prior to the labeled expiration date. Multiple thaw/freeze cycles should be avoided as they will hasten the degradation of serum nutrients and can result in the formation of insoluble precipitates.

United States

SAFC Biosciences
13804 W. 107th Street
Lenexa, Kansas 66215
USA
Phone +1 913-469-5580
Toll free-USA 1 800-255-6032
Fax +1 913-469-5584

Europe

SAFC Biosciences
Smeaton Road, West Portway
Andover, Hampshire SP10 3LF
UNITED KINGDOM
Phone +44 (0)1264-333311
Fax +44 (0)1264-332412

Asia Pacific

SAFC Biosciences
18-20 Export Drive
Brooklyn, Victoria 3025
AUSTRALIA
Phone +61 (0)3-9362-4500
Toll free-AUS 1 800-200-404
Fax +61 (0)3-9315-1656

Preparation Instructions

Thawing

1. Remove the serum bottles from the freezer and allow them to acclimate to room temperature for approximately 10 minutes.
2. Place each container in a 30 to 37 C water bath or incubator. Excessive temperatures will degrade heat labile nutrients. If using a water bath, prevent the bottle caps from being completely submerged.
3. Gently swirl or shake the bottles every 10 - 15 minutes until the serum is completely thawed.
4. After thawing, use the serum promptly. Liquid serum may be stored refrigerated (2 to 8 C) up to four weeks. To avoid thaw/freeze cycles or long periods of refrigeration, it is recommended that any unused serum be immediately dispensed into small, useful aliquots and refrozen for future use.

Periodic agitation is crucial to its optimum performance. If a bottle of serum is not periodically shaken or swirled as it thaws, gradients containing high concentrations of salts, proteins and lipids will form throughout the liquid portion and lead to the formation of crystalline or flocculent precipitates. These cryoprecipitates are not toxic to cell cultures, but they affect the appearance and consistency of each bottle of serum. Small amounts of cryoprecipitates are not uncommon, and will not affect product performance. Gently warming and mixing the serum will generally allow the material to go back into solution.

Characteristics

Adventitious Viral Agents (AVA) (9CFR 113.53)

None detected

Electrophoretic Profile

Normal pattern

Endotoxin

≤ 10.0 EU/mL

Hemoglobin

≤ 20 mg/dL

Mycoplasma (9CFR 113.28)

None detected

Osmolality

260 - 330 mOsm/kg H₂O

pH (at 25 C)

6.8 - 8.1

Sterility (Current USP)

No microbial growth detected

Total Protein

3.0 - 4.3 g/dL

Test results are recorded For Information Only on the following: Chemical Profile, Cloning Efficiency, Growth Promotion, Plating Efficiency and Virus Antibody.

Warranty, Limitation of Remedies

SAFC Biosciences warrants to the purchaser for a period of one year from date of delivery that this product conforms to its specifications. Other terms and conditions of this warranty are contained in SAFC Biosciences' written warranty, a copy of which is available upon request. ALL OTHER WARRANTIES, EXPRESSED OR IMPLIED, INCLUDING THE IMPLIED WARRANTY OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE, ARE EXCLUDED. In no case will SAFC Biosciences be liable for any special, incidental, or consequential damages arising out of this product or the use of this product by the customer or any third party based upon breach of warranty, breach of contract, negligence, strict tort, or any other legal theory. SAFC Biosciences expressly disclaims any warranty against claims by any third party by way of infringement or the like. THIS PRODUCT IS INTENDED FOR PURPOSES DESCRIBED ONLY AND IS NOT INTENDED FOR ANY HUMAN OR THERAPEUTIC USE.

Additional Terms and Conditions are contained in the product Catalog, a copy of which is available upon request.

SER-TAIN™ is a trademark of SAFC Biosciences.

© 2005 SAFC Biosciences

Issued September 2005 XP12106
0805

United States

SAFC Biosciences
13804 W. 107th Street
Lenexa, Kansas 66215
USA
Phone +1 913-469-5580
Toll free-USA 1 800-255-6032
Fax +1 913-469-5584

Europe

SAFC Biosciences
Smeaton Road, West Portway
Andover, Hampshire SP10 3LF
UNITED KINGDOM
Phone +44 (0)1264-333311
Fax +44 (0)1264-332412

Asia Pacific

SAFC Biosciences
18-20 Export Drive
Brooklyn, Victoria 3025
AUSTRALIA
Phone +61 (0)3-9362-4500
Toll free-AUS 1 800-200-404
Fax +61 (0)3-9315-1656