


ChemFiles

Supplement II, 2007



ChiroSolv[®] Resolving Kits

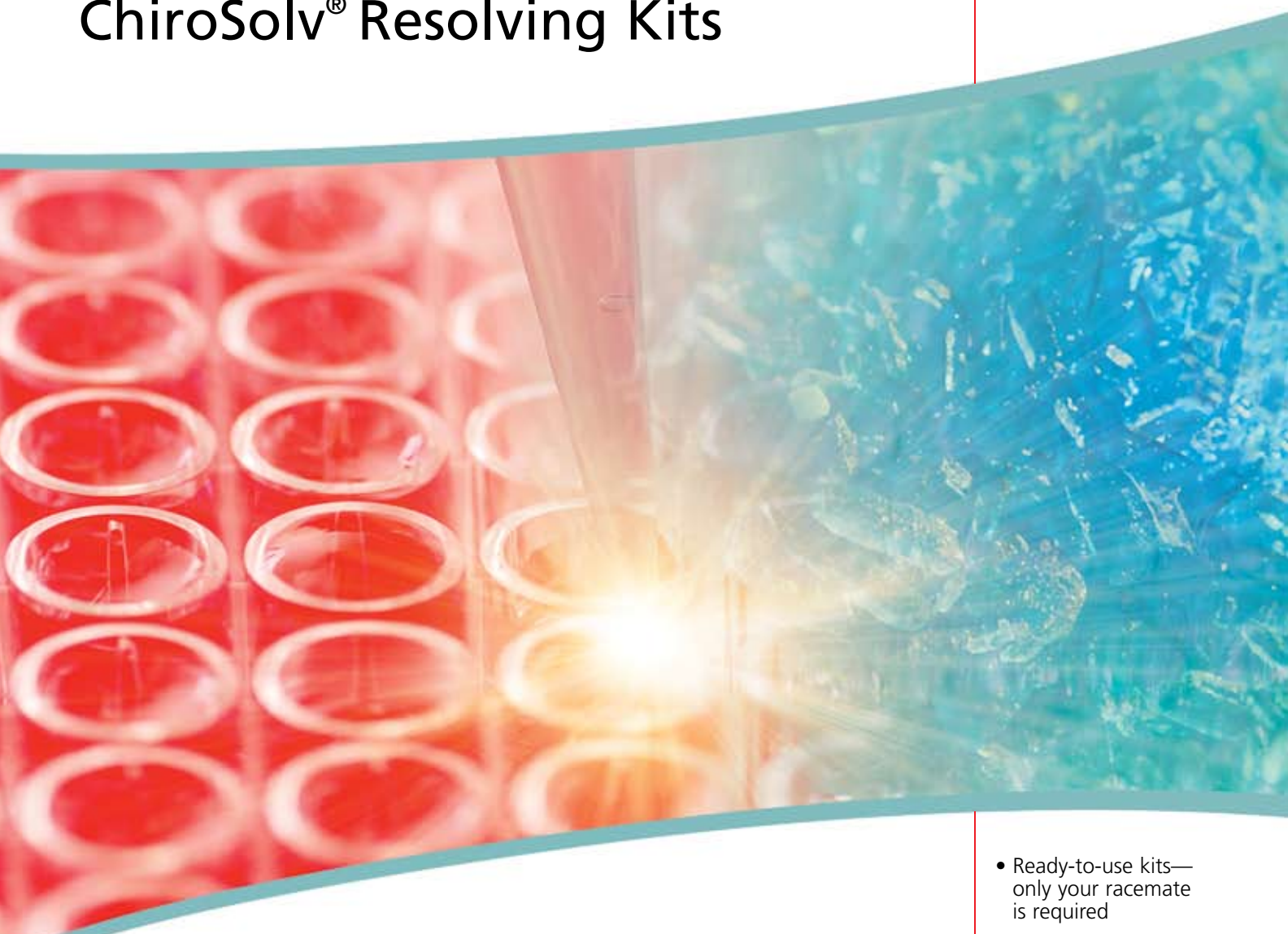

Sigma-Aldrich and ChiroSolve, Inc. are partnering to bring facile and rapid chiral resolution to you.



- Ready-to-use kits—only your racemate is required
- Accurate and consistent results
- Increased likelihood of project success
- Tremendous time and money savings

Facile and Rapid Chiral Resolution with ChiroSolv Kits

Sigma-Aldrich is pleased to partner with ChiroSolve, Inc. to offer, through an exclusive distribution agreement, a series of ready-to-use kits for chiral resolution of racemates. These kits allow scientists to quickly screen resolving agents and solvents against a racemate to find the optimum combination, as well as to optimize reaction conditions in order to separate a racemic mixture into its constituent enantiomers. They can be used for racemic acids, bases, alcohols, amino acids, aldehydes, and ketones.

There are three Acid Series kits and three Base Series kits, available individually or as the complete acid or base set. Each individual kit is in a ready-to-use, 96-vial, high-throughput format that allows the entire analysis to be performed inside the kit without removing any vials. The kit is made of polypropylene material that can withstand extreme temperatures (-20 to 120 °C), allowing the entire kit to be placed in an oven or ice bath without damage.

Each vial has calibrated quantities of a unique resolving agent and solvent combination. Resolving agents are chosen with manufacturing use in mind; they are relatively inexpensive and recoverable in high yield after the resolution is complete. Using all three kits in an acid or base series, 288 resolving agent-solvent combinations can be explored for a single racemate. The high-throughput format allows scientists to identify within one day the optimum resolution process which might otherwise take over 2 months.

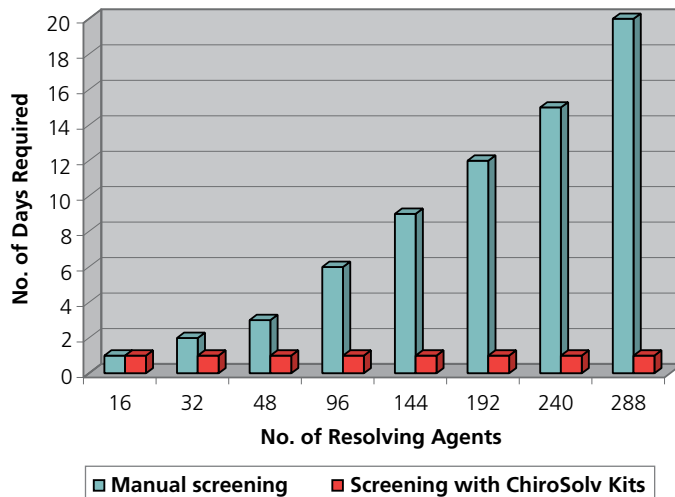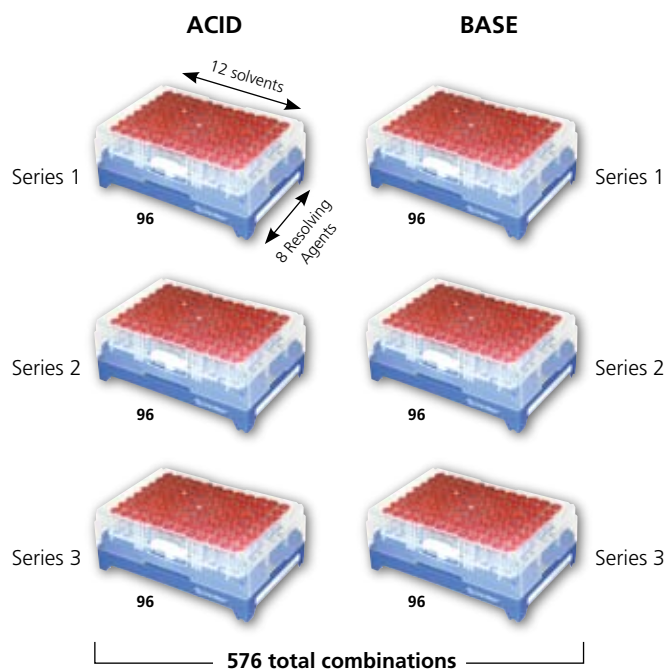

Types of Kits

Acid

Acid Series 1, 2, and 3 include a distinct group of chirally pure acids and are used to resolve racemic bases and amino acids (some pre-processing is required for amino acids). Examples of resolving agents in the acid kits are:

- (-)-Camphoric acid
- (+)- and (-)-Camphorsulphonic acid
- (+)- and (-)-Dibenzoyltartaric acid
- (-)-Malic acid
- (+)- and (-)-Mandelic acid
- (+)-Lactic acid
- (+)- and (-)-Tartaric acid

Base

Base Series 1, 2, and 3 include a distinct group of chirally pure bases or amines and are used to resolve racemic acids, alcohols, aldehydes, and ketones (some pre-processing is required for the latter three). Examples of resolving agents in the base kits are:


- (-)-2-Aminobutanol
- (-)-Brucine
- (-)-Cinchonidine
- (+)-Cinchonine
- (+)-Dehydroabietylamine
- (+)- and (-)-Methylbenzylamine
- (+)-Quinidine
- (-)-Quinine

Case Studies with ChiroSolv Kits


ChiroSolv Kits have been used by several pharmaceutical, fine chemical, and contract R&D companies to reduce their product development time and increase speed to market. Two examples of the application of this technology for determining the ideal solvent/resolving agent combination for the resolution of key compounds are shown.

In the first case, a key advanced intermediate in the synthesis of (*S*)-(-)-cetirizine, a non-sedative histamine H-1 receptor antagonist used for the treatment of allergies, was screened against the kits in the Acid Series. The combination of L-(+)-tartaric acid in methanol was determined to provide the maximum optical purity after its chiral resolution.

In the second example, (-)-ephedrine in isobutanol was determined to be the ideal combination for resolving indoline-2-carboxylic acid to yield the (*S*)-(-)-enantiomer, which is an intermediate for ACE inhibitors such as trandolapril and perindopril. For this compound, the racemic mixture was screened against the three ChiroSolv Kits in the Base Series.

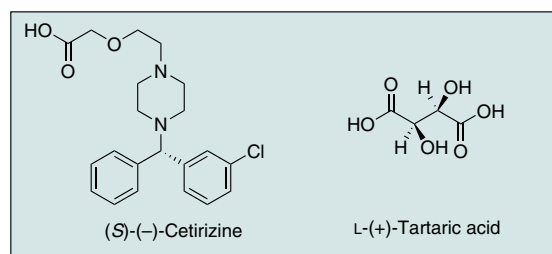

Case Study 1

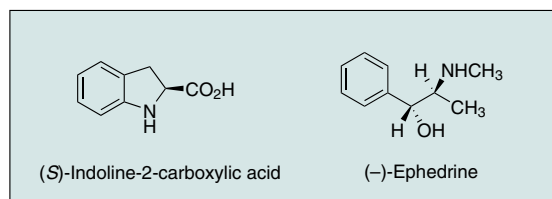

Case Study 2

How Do ChiroSolv Kits Work?

ChiroSolv Kits use the method of diastereomeric crystallization, a process that chemically separates enantiomers in a racemic mixture by complexation with an enantiopure acid or base, resulting in a mixture of diastereomeric salts. These salts have different chemical and physical properties, which allow their separation. This well-established technology competes favorably with newer techniques such as asymmetric synthesis, biocatalysis, enzyme resolution, kinetic resolution, or chiral chromatography.


One major advantage of diastereomeric crystallization is that the procedure scales up easily for manufacturing purposes. Today around 65% of chiral products are made using this technique.

Examples:

- Paroxetine is a selective serotonin reuptake inhibitor developed by Novo Nordisk's subsidiary, Ferrosan, and licensed to GlaxoSmithKline. It is sold as the hydrochloride salt under the name Paxil® for the treatment of depression,

anxiety, and obsessive-compulsive disorder. It is prepared by resolution using (–)-di-*p*-toluoyl-L-tartaric acid.

- Setraline is an antidepressant that inhibits the uptake of serotonin in the central nervous system and is sold by Pfizer under the name Zoloft®. Diastereomeric crystallization of the racemate by (*R*)-mandelic acid is used in its preparation.
- 2-Amino-5-methoxytetralin is an intermediate in the synthesis of N-0923 (Nagase & Co., Ltd., Japan), a potent dopamine D₂ agonist, effective against Parkinson's disease (currently in clinical trial). This intermediate is resolved by diastereomeric salt formation with (*S*)-mandelic acid. The undesired isomer is then racemized and reused in an iterative process.
- In the synthesis of AG-7404 (Pfizer), a key intermediate was resolved by (–)-norephedrine. The final compound was obtained by the combination of diastereomeric and enzymatic resolution.

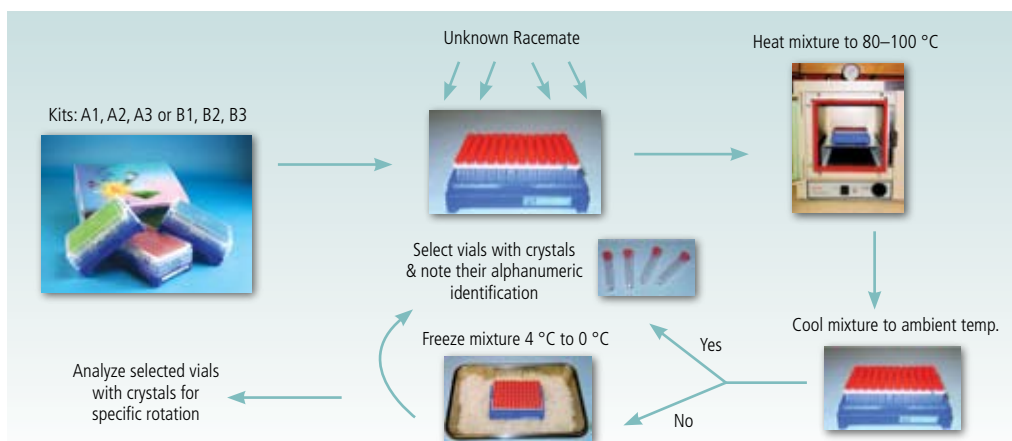


How to Use ChiroSolv Kits

1. Choose the right type of kit (Acid Series 1, 2, 3 or Base Series 1, 2, 3) depending on the identity of the racemate.
2. If the racemate is an alcohol, amino acid, aldehyde or ketone, you may need to pre-process the racemate (refer to <http://www.chirosolve.com/typesofkit.html> for details).
3. Add 0.03 mmol of your racemate to each of the 96 vials. Depending on the availability of the dispensing autostation and the racemate type (liquid or powder), you may need to remove the vial caps. Note that the caps are pierceable to accommodate direct injection of racemate by liquid dispensers.
4. In an oven or water bath, heat the rack along with its vials to 80 °C (optimum temperature for most of these experiments) or until the mixture becomes homogeneous.
5. Allow the kit to cool to ambient temperature. If required, further cool it to 4 °C or 0 °C to observe crystallization. Vials with crystals are considered positive results and should be examined further.
6. Often, more vials with positive results can be obtained using standard crystal initiation techniques. Vials with no crystals even after this effort are considered negative results.
7. Separate out the positive test results and note down their alphanumeric identification.
8. Liberate the enantiomers from their diastereomeric salts and analyze by optical rotation. Typically, only one or two vials will show maximum optical purity.

Features

- Six types of ready-to-use disposable kits providing a total of 576 resolving agent–solvent combinations.
- Very little amount of racemate required (less than 3 mmol per kit).
- Concentrations of the resolving agents are calibrated to give accurate results and optimal crystallization.
- Pierceable rubber septa allow easy racemate dispensing; vials held together by a plastic ring so that vials can be picked up and inspected together for crystal formation.
- Easy to follow, step-by-step instructions and results charts.
- Conveniently designed for easy robotic manipulations.
- Kits and vials have unique identification (barcode or alphanumeric) for cross-referencing.
- Long shelf life at room temperature (2 years).

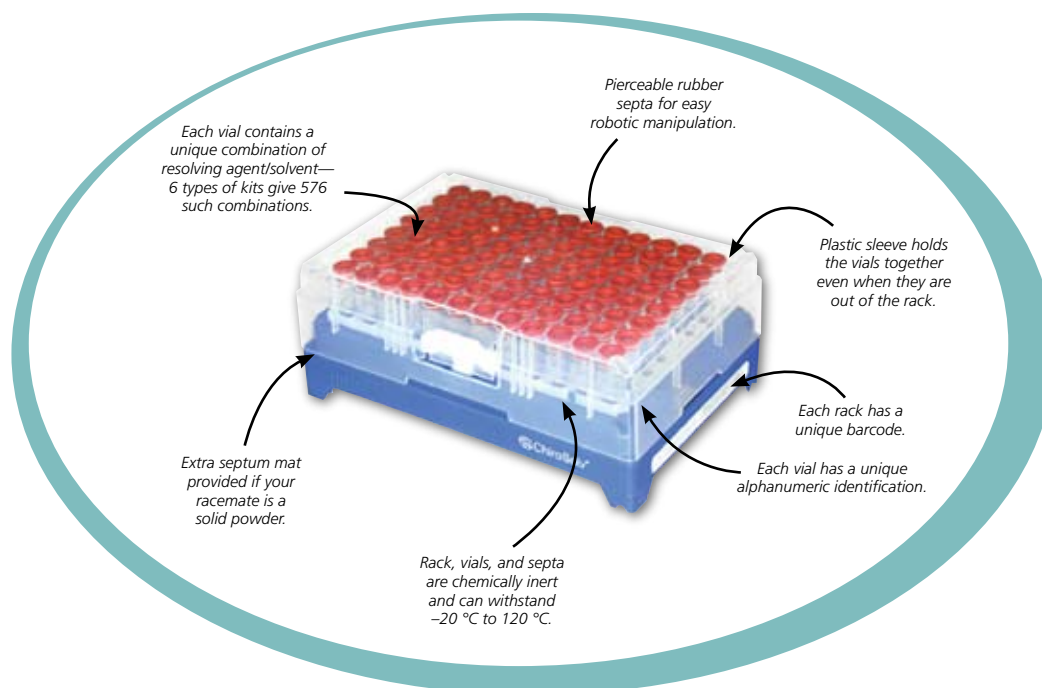

Product Information

Prod. No.	Product Name/Description
681431	ChiroSolv Resolving Kit, Acid Series 1
681423	ChiroSolv Resolving Kit, Acid Series 2
681415	ChiroSolv Resolving Kit, Acid Series 3
688460	ChiroSolv Resolving Kit, Complete Acid Series
681407	ChiroSolv Resolving Kit, Base Series 1
681393	ChiroSolv Resolving Kit, Base Series 2
681377	ChiroSolv Resolving Kit, Base Series 3
688479	ChiroSolv Resolving Kit, Complete Base Series

For more information, please visit sigma-aldrich.com/chirosolv

Argentina

SIGMA-ALDRICH DE ARGENTINA S.A.
Free Tel: 0810 888 7446
Tel: (+54) 11 4556 1472
Fax: (+54) 11 4552 1698

Australia

SIGMA-ALDRICH PTY LTD.
Free Tel: 1800 800 097
Free Fax: 1800 800 096
Tel: (+61) 2 9841 0555
Fax: (+61) 2 9841 0500

Austria

SIGMA-ALDRICH HANDELS GmbH
Tel: (+43) 1 605 81 10
Fax: (+43) 1 605 81 20

Belgium

SIGMA-ALDRICH NV/SA.
Free Tel: 0800 14747
Free Fax: 0800 14745
Tel: (+32) 3 899 13 01
Fax: (+32) 3 899 13 11

Brazil

SIGMA-ALDRICH BRASIL LTDA.
Free Tel: 0800 701 7425
Tel: (+55) 11 3732 3100
Fax: (+55) 11 5522 9895

Canada

SIGMA-ALDRICH CANADA LTD.
Free Tel: 1800 565 1400
Free Fax: 1800 265 3858
Tel: (+1) 905 829 9500
Fax: (+1) 905 829 9292

China

SIGMA-ALDRICH (SHANGHAI)
TRADING CO. LTD.
Free Tel: 800 819 3336
Tel: (+86) 21 6141 5566
Fax: (+86) 21 6141 5567

Czech Republic

SIGMA-ALDRICH spol. s r. o.
Tel: (+420) 246 003 200
Fax: (+420) 246 003 291

Denmark

SIGMA-ALDRICH DENMARK A/S
Tel: (+45) 43 56 59 10
Fax: (+45) 43 56 59 05

Finland

SIGMA-ALDRICH FINLAND OY
Tel: (+358) 9 350 9250
Fax: (+358) 9 350 92555

France

SIGMA-ALDRICH CHIMIE S.à.r.l.
Free Tel: 0800 211 408
Free Fax: 0800 031 052
Tel: (+33) 474 82 28 00
Fax: (+33) 474 95 68 08

Germany

SIGMA-ALDRICH CHEMIE GmbH
Free Tel: 0800 51 55 000
Free Fax: 0800 64 90 000
Tel: (+49) 89 6513 0
Fax: (+49) 89 6513 1160

Greece

SIGMA-ALDRICH (O.M.) LTD.
Tel: (+30) 210 994 8010
Fax: (+30) 210 994 3831

Hungary

SIGMA-ALDRICH Kft
Ingyenes zöld telefon: 06 80 355 355
Ingyenes zöld fax: 06 80 344 344
Tel: (+36) 1 235 9055
Fax: (+36) 1 235 9050

India

SIGMA-ALDRICH CHEMICALS
PRIVATE LIMITED
Telephone
Bangalore: (+91) 80 6621 9600
New Delhi: (+91) 11 4165 4255
Mumbai: (+91) 22 2570 2364
Hyderabad: (+91) 40 4015 5488
Fax
Bangalore: (+91) 80 6621 9650
New Delhi: (+91) 11 4165 4266
Mumbai: (+91) 22 2579 7589
Hyderabad: (+91) 40 4015 5466

Ireland

SIGMA-ALDRICH IRELAND LTD.
Free Tel: 1800 200 888
Free Fax: 1800 600 222
Tel: (+353) 1 404 1900
Fax: (+353) 1 404 1910

Israel

SIGMA-ALDRICH ISRAEL LTD.
Free Tel: 1 800 70 2222
Tel: (+972) 8 948 4100
Fax: (+972) 8 948 4200

Italy

SIGMA-ALDRICH S.r.l.
Numero Verde: 800 827018
Tel: (+39) 02 3341 7310
Fax: (+39) 02 3801 0737

Japan

SIGMA-ALDRICH JAPAN K.K.
Tel: (+81) 3 5796 7300
Fax: (+81) 3 5796 7315

Korea

SIGMA-ALDRICH KOREA
Free Tel: (+82) 80 023 7111
Free Fax: (+82) 80 023 8111
Tel: (+82) 31 329 9000
Fax: (+82) 31 329 9090

Malaysia

SIGMA-ALDRICH (M) SDN. BHD
Tel: (+60) 3 5635 3321
Fax: (+60) 3 5635 4116

Mexico

SIGMA-ALDRICH QUÍMICA, S.A. de C.V.
Free Tel: 01 800 007 5300
Free Fax: 01 800 712 9920
Tel: 52 722 276 1600
Fax: 52 722 276 1601

The Netherlands

SIGMA-ALDRICH CHEMIE BV
Free Tel: 0800 022 9088
Free Fax: 0800 022 9089
Tel: (+31) 78 620 5411
Fax: (+31) 78 620 5421

New Zealand

SIGMA-ALDRICH NEW ZEALAND LTD.
Free Tel: 0800 936 666
Free Fax: 0800 937 777
Tel: (+61) 2 9841 0555
Fax: (+61) 2 9841 0500

Norway

SIGMA-ALDRICH NORWAY AS
Tel: (+47) 23 17 60 60
Fax: (+47) 23 17 60 50

Poland

SIGMA-ALDRICH Sp. z o.o.
Tel: (+48) 61 829 01 00
Fax: (+48) 61 829 01 20

Portugal

SIGMA-ALDRICH QUÍMICA, S.A.
Free Tel: 800 202 180
Free Fax: 800 202 178
Tel: (+351) 21 924 2555
Fax: (+351) 21 924 2610

Russia

SIGMA-ALDRICH RUS, LLC
Tel: +7 (495) 621 6037
+7 (495) 621 5828
Fax: +7 (495) 621 5923

Singapore

SIGMA-ALDRICH PTE. LTD.
Tel: (+65) 6779 1200
Fax: (+65) 6779 1822

South Africa

SIGMA-ALDRICH
SOUTH AFRICA (PTY) LTD.
Free Tel: 0800 1100 75
Free Fax: 0800 1100 79
Tel: (+27) 11 979 1188
Fax: (+27) 11 979 1119

Spain

SIGMA-ALDRICH QUÍMICA, S.A.
Free Tel: 900 101 376
Free Fax: 900 102 028
Tel: (+34) 91 661 99 77
Fax: (+34) 91 661 96 42

Sweden

SIGMA-ALDRICH SWEDEN AB
Tel: (+46) 8 742 4200
Fax: (+46) 8 742 4243

Switzerland

SIGMA-ALDRICH CHEMIE GmbH
Free Tel: 0800 80 00 80
Free Fax: 0800 80 00 81
Tel: (+41) 81 755 2828
Fax: (+41) 81 755 2815

United Kingdom

SIGMA-ALDRICH COMPANY LTD.
Free Tel: 0800 717 181
Free Fax: 0800 378 785
Tel: (+44) 1747 833 000
Fax: (+44) 1747 833 313
SAFC (UK) Free Tel: 01202 712305

United States

SIGMA-ALDRICH
P.O. Box 14508
St. Louis, Missouri 63178
Toll-Free: 800 325 3010
Toll-Free Fax: 800 325 5052
Call Collect: (+1) 314 771 5750
Tel: (+1) 314 771 5765
Fax: (+1) 314 771 5757

Internet

sigma-aldrich.com

World Headquarters

3050 Spruce St., St. Louis, MO 63103
(314) 771-5765
sigma-aldrich.com

Order/Customer Service (800) 325-3010 • Fax (800) 325-5052

Technical Service (800) 325-5832 • sigma-aldrich.com/techservice

Development/Bulk Manufacturing Inquiries SAFC™ (800) 244-1173

©2007 Sigma-Aldrich Co. All rights reserved. SIGMA, SAFC, SAFC™, SIGMA-ALDRICH, ALDRICH, FLUKA, and SUPELCO are trademarks belonging to Sigma-Aldrich Co. and its affiliate Sigma-Aldrich Biotechnology, L.P. Sigma brand products are sold through Sigma-Aldrich, Inc. Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip. Chirosoolv is a registered trademark of Chirosoolve, Inc. Paxil is a registered trademark of GlaxoSmithKline. Zolofit is a registered trademark of Pfizer.

JVV
03568-504740
0097

