

## Product Information

**Monoclonal ANTI-FLAG® M2, Clone M2**  
produced in mouse, purified immunoglobulin

Catalog Number **F3165**

### Product Description

Monoclonal ANTI-FLAG M2 is a purified immunoglobulin, IgG1, monoclonal antibody, isolated from murine ascites fluid, that binds to FLAG® fusion proteins.<sup>1</sup> Unlike ANTI-FLAG M1 antibody, the M2 antibody will recognize the FLAG sequence at the N-terminus, Met-N-terminus, C-terminus, or at an internal site of FLAG fusion proteins. Monoclonal ANTI-FLAG M2 is useful for identification and capture of FLAG fusion proteins by common immunological procedures such as Western blots and immunoprecipitation. It is also useful for affinity purification of FLAG fusion proteins when bound to a solid support.

Monoclonal ANTI-FLAG M2 binding is not calcium dependent.

### Reagent

Supplied in 10 mM sodium phosphate, 150 mM NaCl, pH 7.4, containing 0.02% sodium azide.

### Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

### Preparation Instructions

Dilute the antibody solution from 0.5–10 µg/mL in Tris Buffered Saline, pH 8.0, with 3% nonfat milk, Catalog Number T8793. Adjust the antibody concentration to maximize detection sensitivity and to minimize background.

### Storage/Stability

Store the undiluted antibody at –20 °C in working aliquots. Repeated freezing and thawing is not recommended.

**Note:** Overtime, small amounts of purified antibodies can precipitate from solution due to intermolecular hydrophobic interactions. If a precipitate is observed in this product, briefly centrifuge the vial to pellet the precipitate. Withdraw the desired volume of antibody solution from the clear supernatant for use. This should not alter the performance of the purified antibody in Western blot or immunoprecipitation applications.

### Procedure

Improved Western Blot Method for Detecting FLAG Fusion Proteins using Monoclonal ANTI-FLAG M2.

1. Separate FLAG fusion proteins from sample lysates using a standard sodium dodecyl sulfate polyacrylamide gel electrophoresis (SDS-PAGE) protocol. Load 2.5–10 µg of total lysate protein per lane.
2. Transfer proteins from the gel to an Immobilon®-P or other polyvinylidene difluoride (PVDF) membrane. Nitrocellulose membranes can be used, but typically result in less sensitivity.
3. Wash the blot in at least 0.5 mL/cm<sup>2</sup> of Milli-Q® water for 2–3 minutes with mild agitation.
4. Block the blot with at least 0.5 mL/cm<sup>2</sup> of Tris Buffered Saline, pH 8.0, with 3% nonfat milk, or 50 mM Tris, 0.138 M NaCl, 2.7 mM KCl, pH 8.0, containing 30 mg/mL nonfat dry milk, for 30 minutes at room temperature with agitation (about 50–60 rpm).
5. Remove the blocking agent and wash once with 0.5 mL/cm<sup>2</sup> of TBS, Catalog Number T6664.
6. Add Monoclonal ANTI-FLAG M2 to a final concentration of 10 µg/mL to the blot in at least 0.5 mL/cm<sup>2</sup> of TBS with 3% nonfat dry milk and incubate at room temperature for 30 minutes.  
**Note:** Using less Monoclonal ANTI-FLAG M2 antibody may help to reduce background and cross-reactivity. See the Troubleshooting Guide.

7. Remove the Monoclonal ANTI-FLAG M2 solution and wash once with at least 0.5 mL/cm<sup>2</sup> of TBS.
8. Add Anti-Mouse IgG–Peroxidase, Catalog Number A9044 or equivalent, to at least 0.5 mL/cm<sup>2</sup> of TBS with 3% nonfat dry milk. Use the concentrations listed in Table 1. These concentrations are recommended starting concentrations for the antibodies used in the Western blot. Incubate the blots with shaking at room temperature for 30 minutes.

Table 1.  
Antibody Concentrations

ANTI-FLAG M2 primary antibody	Substrate	Secondary antibody concentration
0.5–10 µg/ml	ECL+™	1:80,000
0.5–10 µg/ml	ECL™	1:10,000

9. Wash the blot eight times for a total of 20 minutes in Tris Buffered Saline, with TWEEN® 20, pH 8.0 (50 mM Tris, 0.138 M NaCl, 2.7 mM KCl, pH 8.0, plus 0.05% TWEEN 20), Catalog Number T9039.
10. Develop the blots with the appropriate substrate for 5 minutes.
11. Expose BioMax™ light film to the blot. Exposure times range from 30 seconds to 10 minutes. It is best to do a quick exposure of 10–30 seconds to determine what exposure time is needed. If the signal is too intense even at the short exposure times let the signal decay from 1–8 hours or longer if necessary and then re-expose the film.

#### Immunofluorescence

Monoclonal ANTI-FLAG M2 may be used in immunofluorescent procedures. A typical concentration for use is 20 µg/mL.<sup>2</sup>

#### **Product Profile**

Protein concentration (E<sub>280</sub>): 2.0 – 5.0 mg/mL

Antigenic binding site:

N-Asp-Tyr-Lys-Asp-Asp-Asp-Lys-C

Specificity: Monoclonal ANTI-FLAG M2 detects a single band of protein on a Western blot from an *E. coli* crude cell lysate.

Sensitivity: Monoclonal ANTI-FLAG M2 detects 2 ng of FLAG-BAP™ fusion protein on a dot blot using chemiluminescent detection.

**Note:** In order to obtain best results, it is recommended that each individual user determine working dilution by titration assay.

#### **References**

1. Brizzard, B.L., *et al.*, Immunoaffinity purification of FLAG epitope-tagged bacterial alkaline phosphatase using a novel monoclonal antibody and peptide elution. *BioTechniques*, **16**, 730-735 (1994)
2. Ciaccia, A.V., and Price, E.M., *IBI FLAG Epitope*, **1**, 4-5 (1992)
3. Bjerrum, O.J., and Heegaard, N.H.H., CRC Handbook of Immunoblotting of Proteins, Volume I, Technical Descriptions, CRC Press, (1988) p. 229-236
4. Dunbar, B.S. (ed.) Protein Blotting: A Practical Approach, IRL Press, NY, p. 67-70 (1994)
5. Fortin, A., *et al.*, A 56- to 54-kilodalton *non grata* signal in immunoblot analysis using the horseradish peroxidase chemiluminescence system. *Biochem. Cell Biol.*, **72**, 239-243 (1994)
6. Harlow, E., and Lane, D., Antibodies: A Laboratory Manual, Cold Spring Harbor Laboratory, Cold Spring Harbor, NY, 1988, Product Code A2926
7. Pampori, N.A., *et al.*, Dilution of the chemiluminescence reagents reduces the background noise on western blots. *BioTechniques*, **18**, 589-590 (1995)

FLAG and ANTI-FLAG are registered trademarks of Sigma-Aldrich® Biotechnology LP and Sigma-Aldrich Co.

The product designations of pFLAG, p3XFLAG, pFLAG-1, pFLAG-2, pFLAGSHIFT, pFLAG-CTS, pFLAG-ATS, pFLAG-MAC, pFLAG-CMV, YEpFLAG, and FLAG-BAP are trademarks of Sigma-Aldrich Biotechnology LP and Sigma-Aldrich Co..

BioMax is a trademark of Eastman Kodak Company  
ECL and ECL+ are trademarks of Amersham Biosciences Ltd.

Immobilon and Milli-Q are registered trademarks of the Millipore Corp.

TWEEN is a registered trademark of Croda International PLC

## Troubleshooting Guide

Problem	Possible Cause	Solution
Fusion protein not detected	Protein not expressed	Verify nucleic acid sequence of FLAG in vector construct. If sequence is present, attempt to optimize expression.
	Target protein poorly represented in sample	Positive controls should always be included. If the positive control works, the sample may not contain the FLAG fusion protein of interest or it may be present at concentrations too low to detect. Immunoprecipitation with Monoclonal ANTI-FLAG M2 Affinity Gel, Catalog Number A2220, may be required for low FLAG fusion protein concentrations. Positive controls available from Sigma: <ul style="list-style-type: none"> <li>Amino-terminal FLAG-BAP Fusion Protein, Catalog Number P7582</li> <li>Carboxy-terminal FLAG-BAP Fusion Protein, Catalog Number P7457</li> <li>Amino-terminal Met-FLAG-BAP Fusion Protein, Catalog Number P5975</li> </ul>
	Defective detection reagents	Run appropriate controls to ensure performance. Use 10 ng/lane of a control FLAG-BAP-fusion protein as a positive control. If no signal is obtained with the control, repeat the procedure using a newer lot of antibody-HRP conjugate and freshly prepared reagents.
	Inadequate exposure time using chemiluminescent system	If no signal is seen, expose for longer times. Sigma recommends trying 30 second to 10 minute exposure times.
	Inappropriate film	Switch to film designated for chemiluminescent detection such as Kodak BioMax Light.
	No target protein present on membrane	Verify transfer by visualizing proteins on the membrane using a Ponceau S solution (Catalog Number P7170). If possible, a positive control should always be run to insure components are functioning. Prestained protein markers (e.g. Catalog Numbers C1992 and C4861) may also be used to verify complete transfer.
	Antigen is covered by blocking reagent due to overblocking	Masking of a signal can occur if the blocking reagent (such as casein or gelatin blocking buffers, Catalog Numbers C7594 or G7663, respectively) is used at too high a concentration. A dilution of 1:1 to 1:3 may be done to decrease the concentration. If the problem persists, use TBS with 3% non-fat dry milk, Catalog Number T8793.
	Antibody concentration not optimal	Determine optimal working dilution for ANTI-FLAG antibody by titration. Consider using more antibody if no signal or weak signal is detected. Also, antibody used at too high a concentration can also cause inhibition of signal especially in chemiluminescent detection systems.
Cross-reactivity	Cellular extract concentration is too high	2.5 to 10 µg per lane of total lysate protein is usually enough to obtain a good signal. Load less cellular extract or serially dilute the cell extract to obtain the optimal signal to noise ratio.
	ANTI-FLAG M2 antibody concentration is too high	Dilute ANTI-FLAG M2 antibody from 0.1 to 0.5 µg/ml. Use TBS with 3% non-fat dry milk as diluent.
	Secondary antibody concentration is too high	Sigma recommends initial dilutions of 1:10,000 for ECL and 1:80,000 for ECL+. Further dilutions may be necessary.
	ANTI-FLAG M2 antibody cross-reacts with naturally occurring FLAG-like epitopes	Increasing the temperature to 37 °C during the blocking, binding and wash steps may reduce cross-reactivity. Lysates from mock-transfected controls (transfected with plasmid without insert DNA) will help distinguish the FLAG-fusion proteins from other cross-reacting proteins.

Sigma brand products are sold through Sigma-Aldrich, Inc.  
Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications.  
Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply.  
Please see reverse side of the invoice or packing slip.