

## 91228 Trichloroacetic acid (TCA)

**CAS Number:** 76-03-9

**Product Description:**

Molecular Formula:	C <sub>2</sub> HCl <sub>3</sub> O <sub>2</sub>
Molecular Weight:	163.4 g/mol
Melting Point:	57-58 °C <sup>1</sup>
Boiling Point:	196-197 °C <sup>1</sup>
pH-Wert:	< 1.0 (0.5 M in H <sub>2</sub> O, 25°C)

Trace elemental analyses have been performed on the BioUltra TCA. BioUltra TCA is for applications which require tight control of elemental content. Trichloroacetic acid (TCA) is a reagent that is used for the precipitation of proteins<sup>2,3</sup> and nucleic acids.<sup>4</sup> TCA is also used as a decalcifier and fixative in microscopy. Addition of TCA to a final concentration of 10% (w/v) will precipitate most proteins from solution. A final concentration of 20% (w/v) may be required for smaller proteins of molecular weight less than 20 kDa. The excess TCA can be removed from protein pellets by washes with buffer.<sup>2</sup> A procedure for the recovery of protein from dilute solutions has been reported.<sup>3</sup> For the precipitation of nucleic acids, a 5% solution of ice cold TCA has been used.<sup>4</sup>

**Solubility / Stability:**

A 100% (w/v) solution of TCA can be made by the addition of 227 ml of water to 500 g of TCA.<sup>5</sup> The resulting clear solution will have a density of approximately 1.45 g/ml. Do not use plastic dishes and containers.

Solution storage temperature 2-8 °C

**References:**

1. The Merck Index, 12th ed., Entry # 9756.
2. Protein Purification Methods: A Practical Approach, Harris, E. L. V., and Angal, S., eds., IRL Press (Oxford, UK: 1989), p. 161.
3. Hwang, B. J., and Chu, G., Trichloroacetic Acid Precipitation by Ultracentrifugation to Concentrate Dilute Protein in Viscous Solution. *Biotechniques*, **20(6)**, 982-984 (1996).
4. Molecular Cloning: A Laboratory Manual, 3rd ed., Sambrook, J. F., et al., Cold Spring Harbor Laboratory Press (Cold Spring Harbor, NY: 2001), pp. A8.25-A8.26.
5. Molecular Cloning: A Laboratory Manual, 3rd ed., Sambrook, J. F., et al., Cold Spring Harbor Laboratory Press (Cold Spring Harbor, NY: 2001), p. A1.29.

**Precautions and Disclaimer**

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

The vibrant M and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates.

Detailed information on trademarks is available via publicly accessible resources.

© 2018 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the US and Canada.

