

09978 Ammonium sulfate

CAS number: 7783-20-2

Product Description:

Molecular Formula: $(\text{NH}_4)_2\text{SO}_4$
Molecular Weight: 132.1 g/mol
pH: 5.0-6.0 (25°C, 1 M H₂O)

This product is soluble in water (435 mg/ml, 3.29 M), yielding a clear, colorless solution. Ammonium sulfate saturation charts at 0 °and 25 °C have been reported.¹³

Store at Room Temperature

Trace elemental analyses have been performed on the Biochemika Ultra ammonium sulfate. The Certificate of Analysis provides lot-specific results. Biochemika Ultra ammonium sulfate is for applications which require tight control of elemental content.

Applications:

Ammonium sulfate is a widely used reagent in molecular biology and chromatography. Applications include the precipitation and fractionation of proteins,^{1,14} purification of antibodies,^{9,10} crystallization of proteins,^{2,3,4,15} protein-nucleic acid complexes,⁵ and of nucleic acids.¹⁷ Ammonium sulfate is also widely used in HPLC of proteins, such as in hydrophobic interaction chromatography.^{6,7,8} It is also used for Sepharose chromatography of tRNAs using reverse salt gradients.¹⁶

The use of ammonium sulfate in long PCR buffer, in PCR lysis solution, and in second-strand cDNA synthesis has been reported.¹¹

A protocol for the HPLC analysis of phospholipids using ammonium sulfate has been reported.¹²

References

1. Englund, S., and Seifter, S., Precipitation Techniques. *Methods Enzymol.*, **182**, 285-300 (1990).
2. McPherson, A., Crystallization of macromolecules: general principles. *Methods Enzymol.*, **114**, 112-120 (1985).
3. Giegé, R., and Ducruix, A., in *Crystallization of Nucleic Acids and Proteins*, Ducruix, A., and Giegé, R., eds., Oxford University Press (Oxford, UK: 1999), pp. 1-16.
4. Tessier, P. M., et al., Self-interaction chromatography: a novel screening method for rational protein crystallization. *Acta Crystallogr. D (Biol. Crystallogr.)*, **58**(Pt 10 Pt 1), 1531-1535 (2002).
5. Yaremchuk, A., et al., Improved crystals of *Thermus thermophilus* prolyl-tRNA synthetase complexed with cognate tRNA obtained by crystallization from precipitate. *Acta Crystallogr. D (Biol. Crystallogr.)*, **56**(Pt 2), 197-199 (2000).
6. Mant, C. T., and Hodges, R. S., in *High-Performance Liquid Chromatography of Peptides and Proteins*, Mant, C. T., and Hodges, R. S., eds., CRC Press (Boca Raton, FL: 1991), pp. 437-450.
7. Hyder, S. M., et al., in *High-Performance Liquid Chromatography of Peptides and Proteins*, Mant, C. T., and Hodges, R. S., eds., CRC Press (Boca Raton, FL: 1991), pp. 451-475.



-
8. Jacob, L. R., in Protein Liquid Chromatography, Kastner, M., ed., Elsevier (Amsterdam: 2000), pp. 235-269.
 9. Kent, U. M., Purification of antibodies using ammonium sulfate fractionation or gel filtration. *Methods Mol. Biol.*, **115**, 11-18 (1999).
 10. Page, M., et al., in Basic Protein and Peptide Protocols, Walker, J. M., ed., Humana Press (Totowa, NJ: 1994), pp. 407-432.
 11. Molecular Cloning: A Laboratory Manual, 3rd ed., Sambrook, J., and Russell, D. W., CSHL Press (Cold Spring Harbor, NY: 2001), pp. 6.22, 8.78, 11.43-11.47.
 12. Guan, Z., et al., Separation and quantitation of phospholipids and their ether analogues by high-performance liquid chromatography. *Anal. Biochem.*, **297(2)**, 137-143 (2001).
 13. Green, A. A., and Hughes, W. L., Protein fractionation on the basis of solubility in aqueous solutions of salts and organic solvents. *Methods Enzymol.*, **1**, 67-90 (1955).
 14. T.D. Pollard, *Meth. Enzymol.* 85, 334 (1982)
 15. G.L. Gilliland, D.R. Davies, *Meth. Enzymol.* 104, 370 (1984)
 16. G.W. Hatfield, *Meth. Enzymol.* 59, 215 (1979)
 17. Dock, A.C., et al., Crystallization of transfer ribonucleic acids *Biochimie* 66, 179-201 (1984)

Precautions and Disclaimer:

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

The vibrant M and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. Detailed information on trademarks is available via publicly accessible resources.
© 2018 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the US and Canada.

