Spectroquant® Data Transfer Software

Operating manual

EN

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DE

Table of contents

1	Introduction	4
1.1	Overview	4
1.2	System requirements	4
1.3	Instruction Manual Overview	4
2	Data transmission	5
2.1	The option window	5
2.2	Store location	6
2.2.1	Availability of the Data Transfer	7
2.3	Operator's Software User Interface.....	8
2.3.1	User Interface Window with Data Transfer Connection.....	8
2.3.2	User Interface Window COM Connection.....	9
2.4	Selecting the Saved Data Format	10
2.4.1	Excel Selection	10
2.4.2	Notepad Selection.....	10
2.5	Transmission of Data	11
3	The menu bar	13
3.1	File	13
3.2	Settings.....	13
3.2.1	Settings Data Transfer	13
3.2.2	Settings COM interface	15
4	Software installation.....	16

1 Introduction

1.1 Overview

The Spectroquant® Data Transfer Software is designed to receive data from the instrument and store it either in an Excel-Spreadsheet or as a .txt-file. Depending on the instrument being used for transmission, an infrared interface module (Spectroquant® Data Transfer) or an RS232 computer interface is required. If the latter is not available, an RS232 to USB adapter can also be used.

In the following table the selectable instrument type number 1 and 2 is related to the corresponding system photometer:

Type	System photometer
1	SQ Move 100
2	SQ Multy

1.2 System requirements

Processor	Pentium® 4/M or similar
RAM	512 MB
Screen resolution	1280 x 1024 pixels
Operating system	Windows® XP, Windows Vista, Windows 7
Excel version *	Excel® 2003 or higher
Disk Space	90 MB

* If an Excel version older than 2003 is being used or Excel software has not been installed, no Excel option will be available when the Software is started (see 2.1).

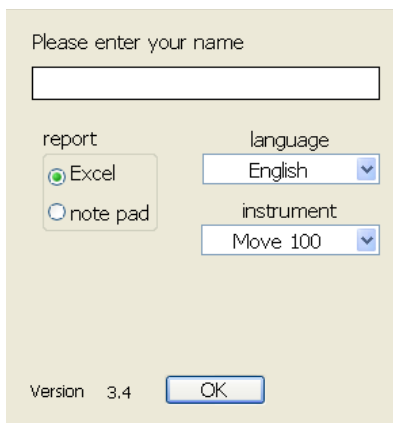
1.3 Instruction Manual Overview

This manual is a reference guide for users of the Data Transfer Software. The user should have prior knowledge of using Windows-based applications. The screenshots provided herein have been created using Windows XP and Excel 2007. If other operating systems or Office versions are being used, differences may occur.

2 Data transmission

2.1 The option window

Once the software has been started, an option window appears. Here you may enter your User Name, select your preferred language and request which type of report is required. You must also select which instrument you are using for the data transmission.

Please enter your name

report

☒ Excel

☐ note pad

language

English


instrument

Move 100

Version 3.4

OK

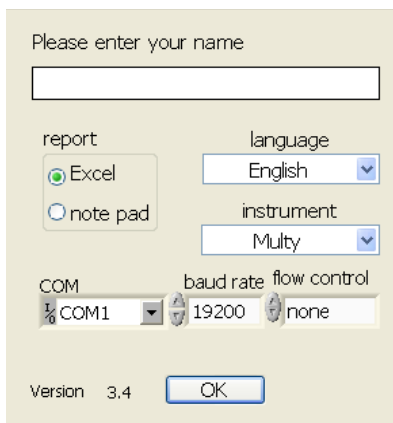If the COM interface is being used for connection, the baud rate and flow control information must be completed. Please also view the instrument's manual for further instructions.

Please enter your name

report

☒ Excel

☐ note pad

language

English

instrument

Multy

COM

COM1

baud rate

19200

flow control


none

Version 3.4

OK

2.2 Store location

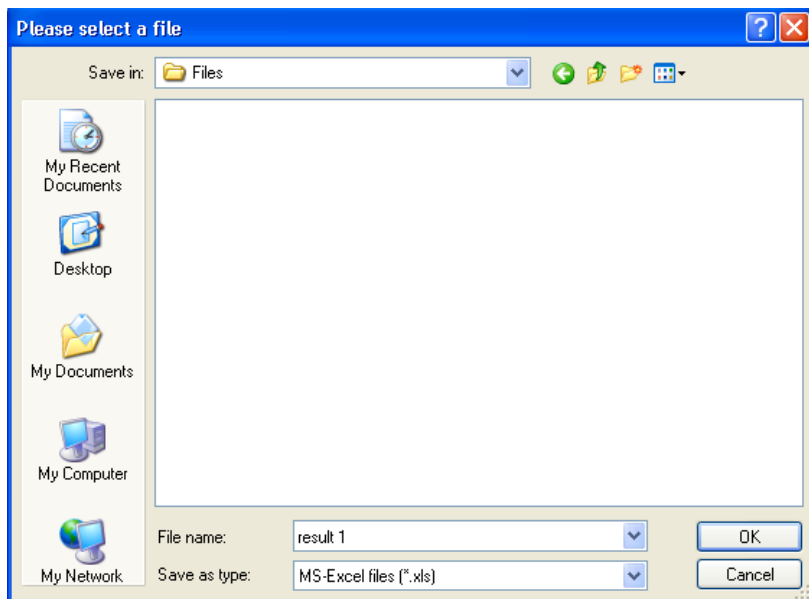
Once the selection has been confirmed, a window appears where the file name and type may be entered. Please note, you may save data either as an Excel Spreadsheet or a Notepad *.txt file.

In Excel, if an existing file name is selected, a new worksheet will be created within this file. If you select to save to an existing Notepad file, the newly transmitted data will be attached to the already existing .txt file.

2.2.1 Availability of the Data Transfer

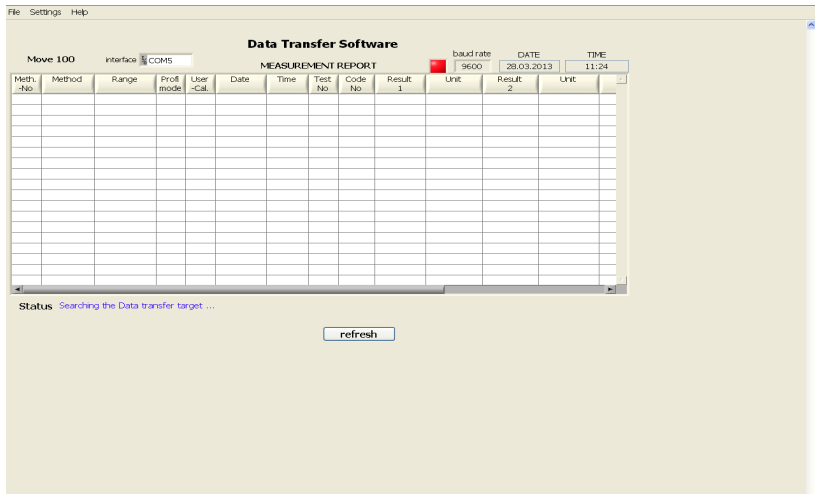After confirming the selection as shown in Section 2.1, the software will automatically verify the Data Transfer's availability. If this verification fails, the following messages will appear in the status box (Example: Move 100).


The screenshot shows the 'Data Transfer Software' window. At the top, there's a menu bar with 'File', 'Settings', and 'Help'. Below it, the 'Move 100' tab is selected. The 'interface' dropdown is set to 'COM5'. The 'MEASUREMENT REPORT' section displays a table with columns: Meth. No., Method, Range, Prof. mode, User Cal., Date, Time, Test No., Code No., Result 1, Unit, Result 2, and Unit. The table is currently empty. To the right of the table, there are fields for 'baud rate' (9600), 'DATE' (28.03.2013), and 'TIME' (11:24). A red status indicator is visible. At the bottom, the 'Status' box shows the message 'Searching the Data transfer target ...' and a 'refresh' button.

Meth. No.	Method	Range	Prof. mode	User Cal.	Date	Time	Test No.	Code No.	Result 1	Unit	Result 2	Unit
-----------	--------	-------	------------	-----------	------	------	----------	----------	----------	------	----------	------

Status: Searching the Data transfer target ...

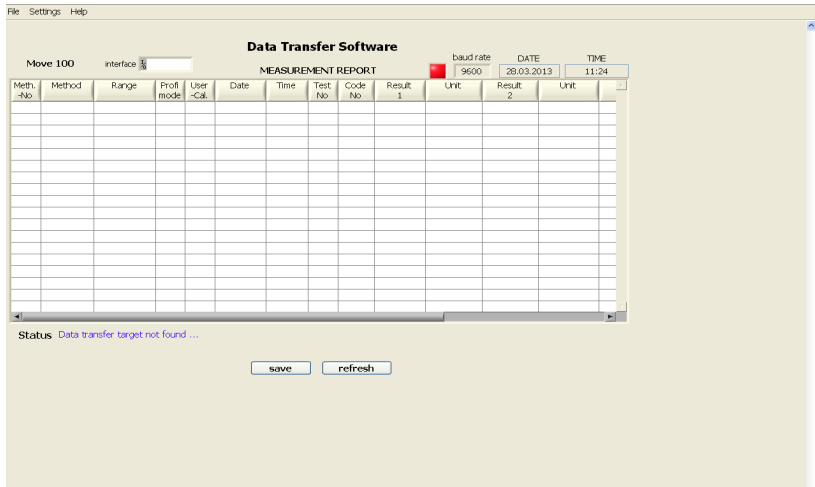refresh

The screenshot shows the 'Data Transfer Software' window with the same layout as the previous one. However, the 'Status' box at the bottom now displays the message 'Data transfer target not found ...'. The 'refresh' button is still present, and a 'save' button has appeared next to it.

Meth. No.	Method	Range	Prof. mode	User Cal.	Date	Time	Test No.	Code No.	Result 1	Unit	Result 2	Unit
-----------	--------	-------	------------	-----------	------	------	----------	----------	----------	------	----------	------

Status: Data transfer target not found ...

save refresh

In this case, the Data Transfer must be reconnected. To establish this has been successful, a new search must be carried out by clicking the Refresh button. Once the Data Transfer has been found to be correctly connected, the storage location window opens. Please confirm the storage location or select an alternative accordingly and press "save".

EN

2.3.1 User Interface Window with Data Transfer Connection

[illegible]

[illegible]

2.4 Selecting the Saved Data Format

The downloaded data can be stored in two kinds of file format: Excel and Notepad as selected under Section 2.1.

2.4.1 Excel Selection

If the Excel format option was selected in Section 2.1, an Excel Template automatically opens when the Save button is pressed. This template will be used to store the transmitted data.

[illegible]

2.4.2 Notepad Selection

If the Notepad format option was selected in Section 2.1, the data will automatically be stored as a .txt file.

2.5 Transmission of Data

The data can now be transmitted from the instrument to the PC. For details on how to do this, please refer to the instrument's Operation Manual. Once transmission has begun, the status row displays "Data receiving". The transmitted data will be displayed in the User Interface Window (Example: Move 100).

File Settings Help

Data Transfer Software

Move 100 interface COM7

MEASUREMENT REPORT

baud rate 9600 DATE 02/04/2013 TIME 08:19

Method No	Method	Range	Prof. mode	User Cal.	Date	Time	Test No	Code No	Result 1	Unit	Result 2
20	Aluminum 14825	20-700 µg/l Al	nen		2013-03-28	11:37:24	1	1	124	µg/l Al	
20	Aluminum 14825	20-700 µg/l Al	nen		2013-03-28	11:38:14	2	2	152	µg/l Al	
111	Calcium 14815	5-160 mg/l Ca	nen		2013-03-28	11:39:40	3	0	59	mg/l Ca	
461	Sulfite 01746	1.0 - 60.0 mg/l SO ₃	nen		2013-03-28	11:40:31	4	0	30.2	mg/l SO ₃	
10	Saurekap. 01758	0.40 - 8.00 mmol/l OH	nen		2013-03-28	11:41:46	5	0	3.64	mmol/l OH	
123	Chlord 01807	0.50 - 5.00 mg/l Cl	nen		2013-03-28	11:43:07	6	1	0.97	mg/l Cl	
123	Chlord 01807	0.50 - 5.00 mg/l Cl	nen		2013-03-28	11:43:59	7	2	1.05	mg/l Cl	

Status Data receiving...

save refresh

On pressing the "save" button, the data will be transmitted to an Excel Spreadsheet where it can be stored.

Data Transfer Software																
Measurement Report										Tested By						
										Device		Move 100				
Method No	Method	Measurement Range	Prof. Mode	User Cal.	Date	Time	Test No	Code No	Result 1	Unit	Result 2	Unit	Result 3	Unit	Result 4	Unit
7	20	Aluminum 14825	20-700 µg/l Al	nen	28.03.13	11:37:24	1	1	124	µg/l Al						
8	20	Aluminum 14825	20-700 µg/l Al	nen	28.03.13	11:38:14	2	2	152	µg/l Al						
9	111	Calcium 14815	5-160 mg/l Ca	nen	28.03.13	11:39:40	3	0	59	mg/l Ca						
10	461	Sulfite 01746	1.0 - 60.0 mg/l SO ₃	nen	28.03.13	11:40:31	4	0	30.2	mg/l SO ₃						
11	10	Saurekap. 01758	0.40 - 8.00 mmol/l OH	nen	28.03.13	11:41:46	5	0	3.64	mmol/l OH						
12	123	Chlord 01807	0.50 - 5.00 mg/l Cl	nen	28.03.13	11:43:07	6	1	0.97	mg/l Cl						
13	123	Chlord 01807	0.50 - 5.00 mg/l Cl	nen	28.03.13	11:43:59	7	2	1.05	mg/l Cl						
14	30	Ammonium 14739	10-2000 µg/l NH ₄ -N	nen	28.03.13	13:35:09	8	5	1764	µg/l NH ₄ -N						

On pressing the "save" button, the data will be transmitted to an Excel In the case of Notepad, the data is automatically stored as a .txt-file.

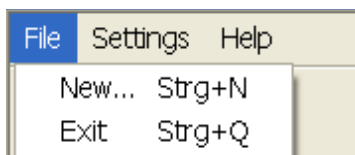
```

Datei Bearbeiten Format Ansicht ?
This record is created by Data Transfer Software
Tested by:
Instrument: Move 100
Record creation date: 28.03.2013
Record creation time: 13:37
-----
20 Aluminium 14825
20-700 µg/l Al
Profil-mode: nein
2013-03-28 11:37:24
laufende Nr.: 1
Code-Nr.: 1
124 µg/l Al

20 Aluminium 14825
20-700 µg/l Al
Profil-mode: nein
2013-03-28 11:38:14
laufende Nr.: 2
Code-Nr.: 2
132 µg/l Al

111 Calcium 14815
5-160 mg/l Ca
Profil-mode: nein
2013-03-28 11:39:40
laufende Nr.: 3
Code-Nr.: 0
59 mg/l Ca

461 Sulfid 01740
1.0 - 50.0 mg/l SO3
Profil-mode: nein
2013-03-28 11:40:31
laufende Nr.: 4
Code-Nr.: 0
30.2 mg/l SO3

```

3 The menu bar

3.1 File


Under the File Tab, the option "New" enables the user to start a new Data Transmission procedure. The software starts again with the selection window shown at 2.1

EN

The software can be closed by using the option File → Exit.

3.2 Settings

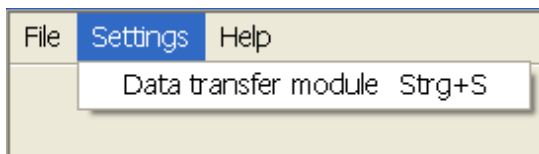
3.2.1 Settings Data Transfer

Under Settings → Data Transfer Module, the Data Transfer's settings can be changed:

1. Baud rate
2. Flow control
3. Automatic switch-off time (only active if the Data Transfer is running on batteries)

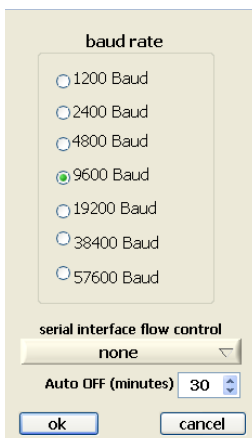
Please see the instrument's manual for further instructions.

The screenshot shows a settings dialog box for the Data Transfer Module. It has a title bar and a main area with the following elements:

- baud rate**: A group box containing seven radio buttons for different baud rates: 1200 Baud, 2400 Baud, 4800 Baud, 9600 Baud (which is selected), 19200 Baud, 38400 Baud, and 57600 Baud.
- serial interface flow control**: A dropdown menu currently set to "none".
- Auto OFF (minutes)**: A numeric input field with a spinner, currently set to 30.
- Buttons**: "ok" and "cancel" buttons at the bottom.

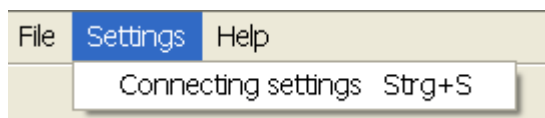
3.2.2 Settings COM interface

The settings of the COM interface can be modified under the Settings Menu → Connecting settings (see also Section 2.1.1):

1. Modify the COM interface to be used
2. Baud rate
3. Flow control

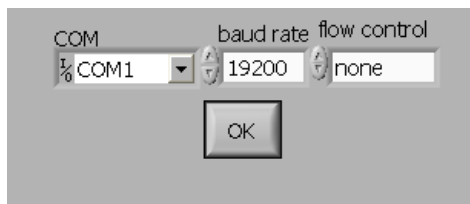
Please see the instrument's manual for further instructions.

4 Software installation


Note: If an older version of the Data Transfer Software is already installed, please remove the program before installing the new software. Please remove National Instrument Software and Data Transfer Software.

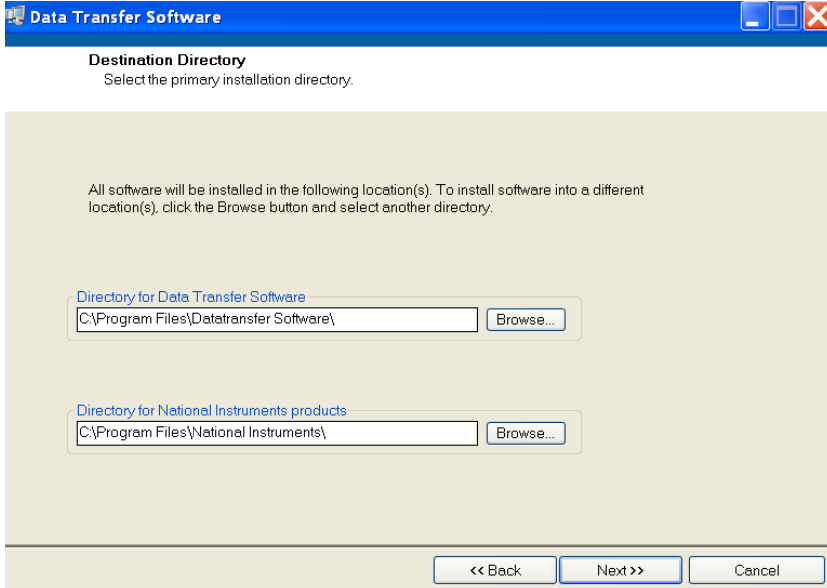On automatic start-up of the setup.exe, the location for program installation has to be confirmed. It is recommended to confirm the default file path by pressing the "Next" button.

On confirmation, two license agreements will appear which must be accepted and confirmed by pressing the "Next" Button.

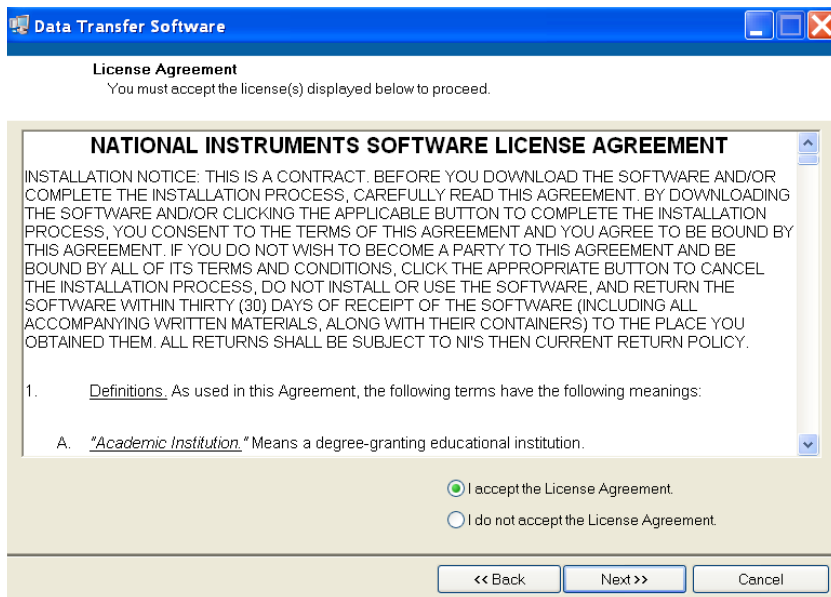On confirmation, two license agreements will appear which must be accepted and confirmed by pressing the "Next" Button.

Data Transfer Software

License Agreement
You must accept the license(s) displayed below to proceed.

NATIONAL INSTRUMENTS SOFTWARE LICENSE AGREEMENT

INSTALLATION NOTICE: THIS IS A CONTRACT. BEFORE YOU DOWNLOAD THE SOFTWARE AND/OR COMPLETE THE INSTALLATION PROCESS, CAREFULLY READ THIS AGREEMENT. BY DOWNLOADING THE SOFTWARE AND/OR CLICKING THE APPLICABLE BUTTON TO COMPLETE THE INSTALLATION PROCESS, YOU CONSENT TO THE TERMS OF THIS AGREEMENT AND YOU AGREE TO BE BOUND BY THIS AGREEMENT. IF YOU DO NOT WISH TO BECOME A PARTY TO THIS AGREEMENT AND BE BOUND BY ALL OF ITS TERMS AND CONDITIONS, CLICK THE APPROPRIATE BUTTON TO CANCEL THE INSTALLATION PROCESS, DO NOT INSTALL OR USE THE SOFTWARE, AND RETURN THE SOFTWARE WITHIN THIRTY (30) DAYS OF RECEIPT OF THE SOFTWARE (INCLUDING ALL ACCOMPANYING WRITTEN MATERIALS, ALONG WITH THEIR CONTAINERS) TO THE PLACE YOU OBTAINED THEM. ALL RETURNS SHALL BE SUBJECT TO NI'S THEN CURRENT RETURN POLICY.

1. Definitions. As used in this Agreement, the following terms have the following meanings:

A. "Academic Institution." Means a degree-granting educational institution.

☒ I accept the License Agreement.
☐ I do not accept the License Agreement.

<< Back Next >> Cancel

Data Transfer Software

License Agreement
You must accept the license(s) displayed below to proceed.

LICENSE AGREEMENT

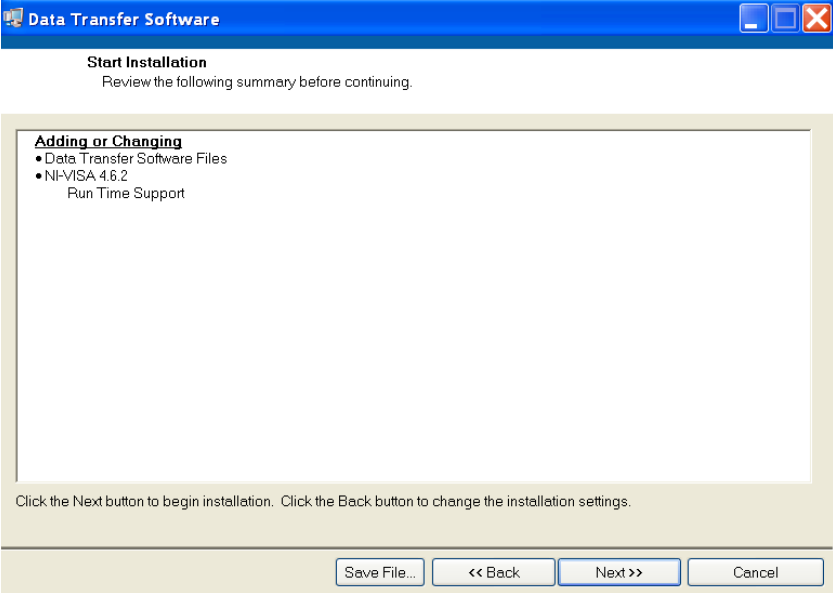
BEFORE YOU CLICK ON THE ACCEPT BUTTON AT THE END OF THIS DOCUMENT, CAREFULLY READ ALL THE TERMS AND CONDITIONS OF THIS AGREEMENT. BY CLICKING ON THE ACCEPT BUTTON, YOU ARE CONSENTING TO BE BOUND BY AND ARE BECOMING A PARTY TO THIS AGREEMENT. IF YOU DO NOT AGREE TO ALL OF THE TERMS OF THIS AGREEMENT, CLICK THE "DO NOT ACCEPT" BUTTON AND DO NOT DOWNLOAD AND/OR USE THIS INTELLECTUAL PROPERTY.

Readers of this document are requested to submit to Interchangeable Virtual Instruments, Inc. ("Licensor"), with their comments, notification of any relevant patent rights or other intellectual property rights of which they may be aware which might be infringed by any use of this intellectual property, software, or specification (the "Intellectual Property"), as appropriate, and to provide supporting documentation.


☒ I accept the License Agreement.
☐ I do not accept the License Agreement.

<< Back Next >> Cancel


Finally, a summary of the program to be installed is displayed. To start the installation press the "Next" button.


The installation starts.

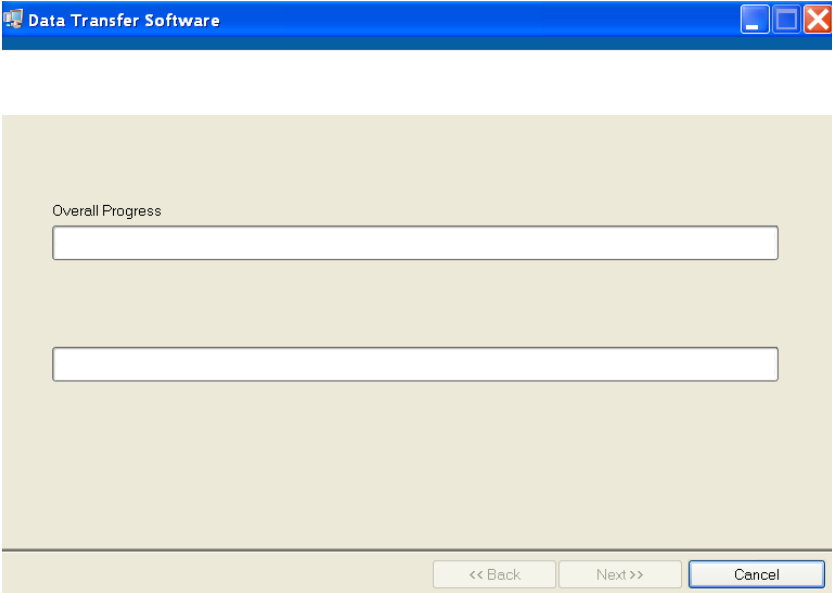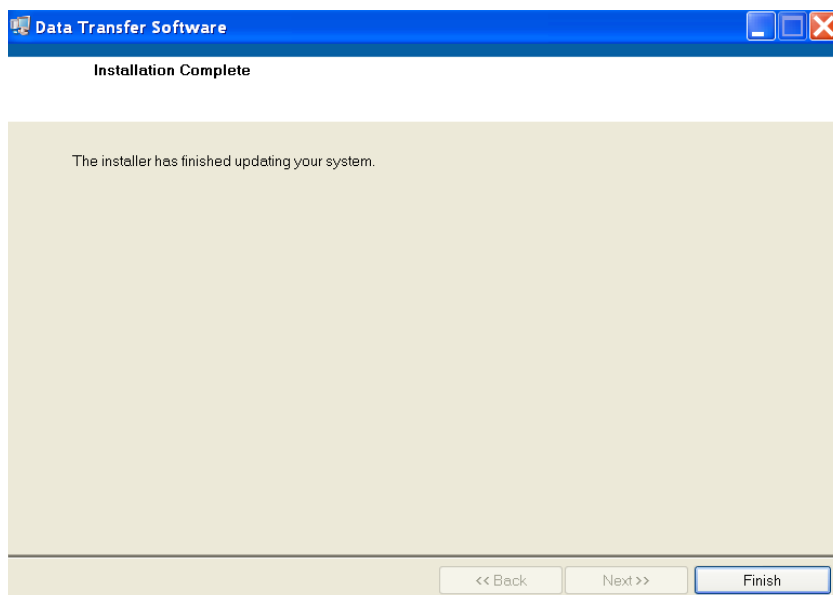Once the installation has been successfully carried out, pressing the "Finish" button ends the procedure.

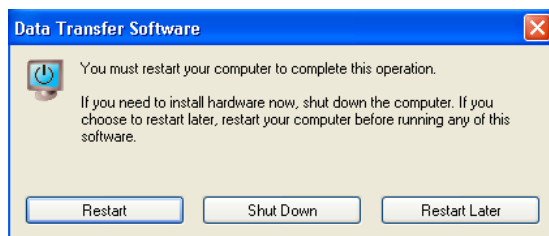After the installation, a Restart of the PC is necessary. On restarting the PC, a link will be created on the desktop which can, in future, be used to start the software.

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
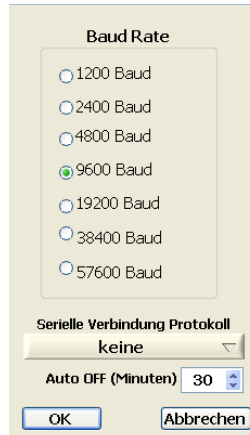
The image shows a settings window titled "Baud Rate". It contains a list of radio buttons for different baud rates: 1200 Baud, 2400 Baud, 4800 Baud, 9600 Baud (which is selected with a green dot), 19200 Baud, 38400 Baud, and 57600 Baud. Below this list is a section titled "Serielle Verbindung Protokoll" with a dropdown menu currently showing "keine". Underneath is a label "Auto OFF (Minuten)" followed by a numeric input field containing "30" and a small up/down arrow. At the bottom of the window are two buttons: "OK" and "Abbrechen".

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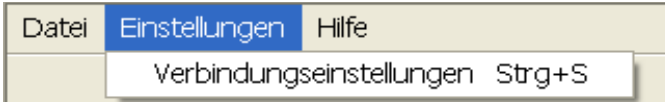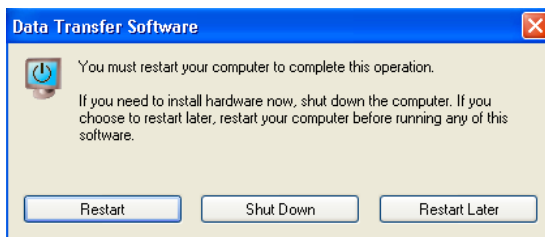
We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

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The vibrant M, Supelco and Spectroquant are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2019 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

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Vertrieb: EMD Millipore Corporation, 400 Summit Drive, Burlington MA 01803, USA
Sigma-Aldrich Canada Co. or Millipore (Canada) Ltd. , 2149 Winston Park,