

Product Information

Apyrase from potato

Catalog Number **A6535**
Storage Temperature $-20\text{ }^{\circ}\text{C}$

CAS RN 9000-95-7
EC 3.6.1.5

Synonyms: Adenosine 5'-diphosphatase, Adenosine 5'-triphosphatase, ATP diphosphohydrolase¹

Product Description

Apyrase has adenosine 5'-triphosphatase (ATPase) and adenosine 5'-diphosphatase (ADPase) activities,² where hydrolysis of the pyrophosphate bonds leads to sequential release of inorganic orthophosphate.³ At least two isoenzymes with different ATPase/ADPase ratios exist in different varieties of potato (*Solanum tuberosum*).^{4,5}

- 'Pimpernel' isoenzyme, with a high ATPase/ADPase ratio of ~10:1 and an isoelectric point (pI) of 8.74
- 'Desirée' isoenzyme, with a low ATPase/ADPase ratio of ~1:1 and an isoelectric point (pI) of 6.69

These isozymes each have a molecular mass of ~49 kDa (gel filtration).⁵ This product is predominantly the low ATPase/ADPase ratio isozyme.

Apyrase requires divalent metal ions for activity, with optimal activity observed with 5 mM Ca^{2+} . For hydrolysis of organic di- and triphosphates, the optimal pH is 6.⁵ For inorganic substrates, the optimal pH is 5.1.⁵

Preparation Instructions

This product is soluble in water (1 mg/mL). One publication reports preparation of 125 units/mL stock solutions of this product in 10 mM HEPES, pH 7.8.⁶

Storage/Stability

Stock solutions at pH between 5–7.5 can be stored as frozen aliquots, such as in 30 mM HEPES, pH 7.2.⁷ For enzyme solutions of <1 mg/ml, dissolve in HEPES buffer, pH 7.5, containing 1 mM MgCl_2 , 1 mM DTT, 1 mM EDTA, and 1 mg/mL BSA. Repeated freeze-thaw cycles risk loss of activity. One publication indicates storage of 340 units/mL stock solutions of apyrase in calcium-free Tyrode's buffer, at $-20\text{ }^{\circ}\text{C}$, in single-use aliquots.⁸

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

References

1. Vasconcelos, E.G., *et al.*, "Potato apyrase: a new tool for parasitic disease research", in *Potato II. Fruit, Vegetable and Cereal Science and Biotechnology 3 - Special Issue 1* (P. Tennant and N. Benkeblia, eds.). Global Science Books, pp. 80-88 (2009).
2. Traverso-Cori, A., *et al.*, *Arch. Biochem. Biophys.*, **109(1)**, 173-184 (1965).
3. Mancilla, M., *et al.*, *Phytochemistry*, **23(7)**, 1397-1400 (1984).
4. Molnar, J., and Lorand, L., *Arch. Biochem. Biophys.*, **93(2)**, 353-363 (1961).
5. Kettlun, A., *et al.*, *Phytochemistry*, **21(3)**, 551-558 (1982).
6. Gonda, K., and Kikyo, N., "Nuclear Remodeling Assay in *Xenopus* Egg Extract", in *Nuclear Transfer Protocols: Cell Programming and Transgenesis* (P.J. Verma and A. Trounson, eds.), *Methods in Molecular Biology*, Vol. 348. Humana Press (Totowa, NJ), Chapter 17, pp. 247-258 (2006).
7. Gennerich, A., and Reck-Peterson, S.L., "Probing the Force Generation and Stepping Behavior of Cytoplasmic Dynein", in *Single Molecule Analysis: Methods and Protocols, Methods in Molecular Biology*, Vol. 783, pp. 63-80 (2011).
8. Rutledge, T.W., and Whitehart, S.W., "Studies of Secretion from Permeabilized Platelets", in *Methods in Molecular Biology: Platelets and Megakaryocytes, Vol. 1 - Functional Assays*, Vol. 272 (J.M. Gibbins and M.P. Mahaut-Smith, eds.). Humana Press (Totowa, NJ), pp. 109-120 (2004).

GCY,MES,NSB,MAM 01/17-1