

## Product Information

# 5-Bromo-4-chloro-3-indolyl $\beta$ -D-glucuronide cyclohexylammonium salt

Tablet

**B8049**

## Product Description

Synonyms: X-GlcA; BC-Indicator; X-glucuronide (5-Bromo-4-chloro-3-indolyl  $\beta$ -D-glucuronide component)

CAS Registry Number: 114162-64-0  
(X-GlcA component)

Molecular Formula:  $C_{14}H_{12}BrClNO_7 \cdot C_6H_{13}N$   
(X-GlcA component)


Molecular Weight: 521.79 (X-GlcA component)

The *gus* operon in *Escherichia coli* is composed of three genes:<sup>1</sup>

- *uidA* (*gusA*), which encodes  $\beta$ -glucuronidase (GUS).
- *gusB*, which encodes a glucuronide permease
- *gusC*, which encodes a 44 kDa protein located in the outer membrane of *E. coli*, with an as-yet undetermined function<sup>2</sup>

5-bromo-4-chloro-3-indolyl  $\beta$ -D-glucuronide (X-GlcA, X-Gluc) has been shown to be a good substrate for GUS, yielding a dark-blue insoluble cleavage product. The reaction (see Figure 1) initially yields a monomeric intermediate, which rapidly oxidizes to form the dimer dichloro-dibromoindigo (Cl Br-indigo).

**Figure 1.** Hydrolysis of X-GlcA by  $\beta$ -Glucuronidase

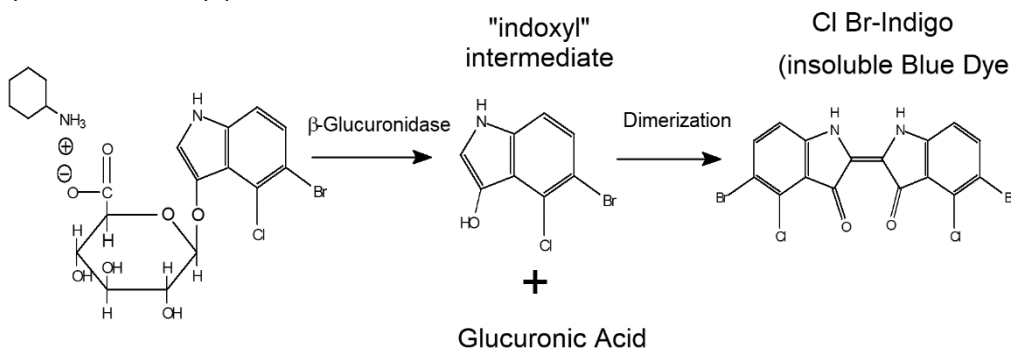

The intense coloration and insolubility of Cl Br-indigo is ideal for use as an indicator of GUS activity *in situ*. Cl Br-indigo has been used as an indicator of *E. coli* contamination in various food items<sup>3</sup> and as an agent in urinary tract infections.<sup>4</sup> The *gusA* gene has been used as an indicator of transfection and as a reporter gene for the function of regulatory elements in plants.<sup>5,6</sup>

If using a known strain of *E. coli* as a positive control for GUS activity, it is important to realize that K-12 strains of *E. coli* contain a defective permease.<sup>1</sup> Even though X-GlcA is an excellent inducer of *uidA* in *E. coli*, K-12 strains require much higher levels of X-GlcA than wild-type strains. With a defective permease, high extracellular levels of X-GlcA are needed to develop sufficient intracellular levels so that *uidA* is adequately induced. In addition, once *uidA* is induced and GUS activity is high, high extracellular levels of X-GlcA are also needed to develop sufficient intracellular levels to react and to yield a dark coloration.

## Reagent

Each tablet is ~40 mg and contains 10 mg of X-GlcA substrate.

## Precautions and Disclaimer

This product is for R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

## Storage/Stability

Store the tablets at –20 °C. When stored at –20 °C, the tablets are stable for at least one year. Tablets are good as long as their color remains white.

## Preparation Instructions

To prepare a concentrated stock solution, dissolve one tablet either in:

- 100 µL of *N,N*-dimethylformamide (DMF) with a final volume of ~125 µL (80 mg/mL final concentration)
- 180 µL of dimethyl sulfoxide (DMSO) with a final volume of 200 µL (50 mg/mL final concentration)

These stock solutions can be frozen at either –20 °C or –70 °C.

## Procedure

As an indicator for the presence of *E. coli* in natural materials, one functional test procedure is as follows:

1. Prepare LB Agar (Cat. No. L2897) or LB Agar EZMix™ Powder (Cat. No. L7533).
2. Cool to 55 °C.
3. Add 250 µL of a 40 mg/mL stock solution of Cat. No. B8049 in DMSO to 100 mL of LB agar. Mix gently to dissolve. The final concentration of X-GlcA in the medium will be 100 µg/mL.
4. Pour plates. Allow to cool for a few hours or overnight.
5. Streak one plate with a *uidA*<sup>+</sup> strain of *E. coli* (ATCC 11303) and a second plate with a *uidA*<sup>-</sup> strain of *E. coli* (GMS407).
6. Incubate the plates at 37 °C for 24 hours.

As a substrate for the GUS reporter system to study plant gene expression, various published procedures are available.<sup>1</sup>

## References

1. Gallagher, S.R. (ed.), *GUS Protocols: Using the GUS Gene as a Reporter of Gene Expression*. Academic Press, Inc. (San Diego, CA: 1992).
2. Liang, W.-J. et al., *J. Bacteriol.*, **187(7)**, 2377-2385 (2005).
3. Delisle, G.J., and Ley, A., *J. Clin. Microbiol.*, **27(4)**, 778-779 (1989).
4. Restaino, L. et al., *J. Food. Prot.*, **53(6)**, 508-510 (1990).
5. Bomineni, V.R. et al., *Plant Cell Rep.*, **13(1)**, 17-23 (1993).
6. Ellis, D.D. et al., *Bio/Technology*, **11**, 84-89 (1993).

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

## Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

## Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

## Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2007-2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

B8049dat Rev 02/23 RBG,GCY,MAM

**MILLIPORE  
SIGMA**