

## Product Information Sheet

# Collagenase from Clostridium histolyticum

Type XI

**C7657**

## Product Description

CAS Number: 9001-12-1

Molecular Weight: This collagenase obtained from the culture filtrate of Clostridium histolyticum. The culture filtrate that is obtained is thought to contain at least 7 different proteases ranging in molecular weight from 68-130 kDa.

Many references exist for using collagenase to digest various tissue. The choice of one technique over another is mostly arbitrary and based more on past experience than on an understanding of why the method works and what modifications could lead to even better results. Concentrations typically vary from 0.1 mg/mL to 5 mg/mL, and digestion time is something that should be experimentally monitored using a very gentle agitation system to check for tissue dissociation. The concentration of collagenase needed to isolate islets is dependent on the method used. A general recommendation is to start at a collagenase concentration of 2.5 mg/mL. Satisfactory dissociation versus viability typically is achieved from 15 minutes to several hours but can fall outside of this range if the concentration is unusual.

The preferred buffer to use is Krebs Ringer Buffer with calcium and BSA. Zn<sup>2+</sup> is required for activity, but it is tightly bound to the collagenase during purification. Additional Zn<sup>2+</sup> should not be necessary as long as no chelator is added to the solution during digestion.

If excessive cell death is observed using concentrations tested in the past, the particular lot used might have a higher specific activity. Lowering the enzyme concentration and/or adding BSA or serum (0.5% and 5-10%, respectively) is recommended. These components are added to stabilize the cells to further digestion by the enzyme.

To sterile filter solutions of collagenase, first centrifuge the solution or filter through a 0.8 µm filter to remove insolubles. This will remove particulates and reduce the probability of clogging the 0.2 µm filter during sterile filtration.

Radiolabeled gelatin has been used to measure the activity and mechanism of collagenase digestion.

Mandl units have the same description as Sigma collagen digestion units. The conversion factor for Mandl units/Wuensch units to Sigma units is approximately 1000-2000 to 1.

Most of the proteases in this preparation may be inhibited by a combination of benzamidine hydrochloride (Cat. No. B6506) and TLCK (Cat. No. T7254) without affecting the activity of collagenase. However, there is no guarantee that all the proteases will be inhibited. For general applications, the working concentrations for benzamidine hydrochloride is about 1 mM and TLCK is 0.100-0.135 mM. If a collagenase with no significant protease activity is required, Cat. No. C0773 is recommended, since this product has been tested for any residual protease activity and found to contain less than 1-unit neutral protease.

## Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

## Preparation Instructions

For in enzymatic activity measurement, an enzyme stock solution is prepared by dissolving 0.05-0.1 mg/mL collagenase in 50 mM TES buffer, pH 7.4, containing 0.36 mM calcium chloride at 37 °C. Final concentrations in the reaction mixture are 50 mM TES, 0.36 mM calcium chloride, 25 mg collagen (Cat. No. C9879), and 0.005-0.01 mg collagenase.

## Storage and Stability

Solutions at neutral pH and with adequate calcium (0.3-0.5 mM) will retain activity for at least 5 hours at 37 °C. Solutions at -20 °C are stable for several months.

## References

1. Angleton, E.L., et. Al., Preparation and reconstitution with divalent metal ions of class I and class II Clostridium histolyticum apocollagenases. *Biochemistry*, 27(19), 7406-7412 (1988).
2. Mookhtiar, K.A., et al., Properties of radiolabeled type I, II, and III collagens related to their use as substrates in collagenase assays. *Anal. Biochem.*, 158(2), 322-333 (1986).
3. Schomberg, et al., *Enzyme Handbook* (Springer-Verlag, Berlin Heidelberg: 1991), p. 1-6

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

### Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

### Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

### Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

C7657pis Rev 08/22

**MILLIPORE  
SIGMA**